


Short-form operating instructions

METTLER TOLEDO SM balances

METTLER TOLEDO

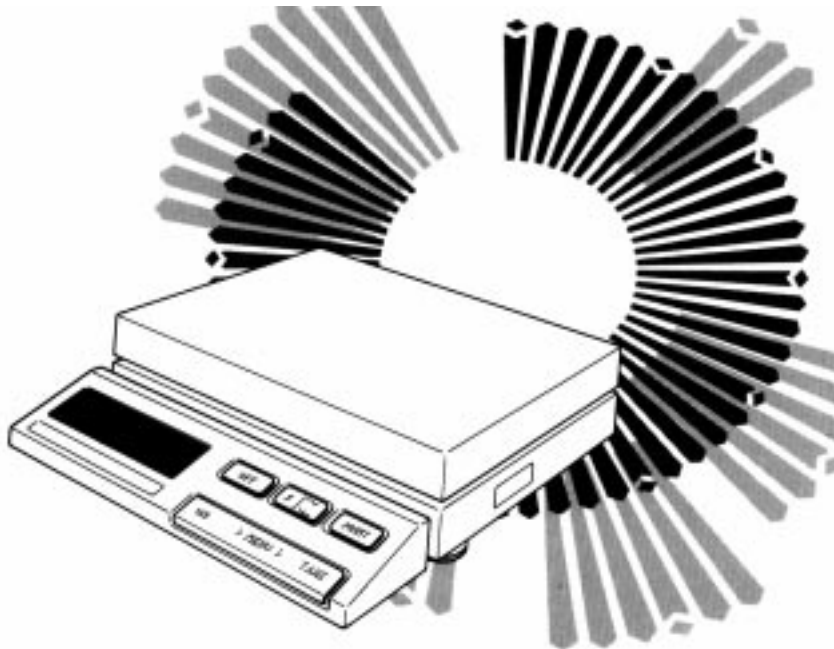

Subject to technical changes.



P703883

© Mettler-Toledo GmbH 1998 703883 Printed in Switzerland 9811/7.12

Mettler-Toledo GmbH, Laboratory & Weighing Technologies CH-8606 Greifensee, Switzerland
Phone +41-1-944 22 11, Fax +41-1-944 30 60, Internet: <http://www.mt.com>


Switching the equipment on and off



Switching on: Briefly press control bar (3).
Note: The scale first goes through an automatic checking sequence and displays 0 <Zero> when ready.
Switching off: Press OFF key (4).

Net / Gross - Weighing


Net

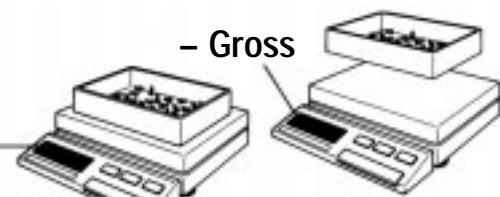

Place empty container on scale and tare.
Fill in → read off net weight.
Tare, then remove everything.
Display → read off - gross weight.


In standard configuration: Plus / minus - percent display




Place weighing goods on pan.
Press F key (5). The display now changes to 100% with the Plus/minus display:
Tolerance limit $\pm 2.5\%$; 100% = nominal weight (minimum weight = 10.0 g)
Plus/minus weighings can now be performed.
Note:
By pressing  key (5) the weight can be read in weight units. Repeated pressing of F key (5) records the weight on the scale as 100%.
The same function key (5) permits still other applications. Please consult the Detailed Operation Instructions.

PRINT Print Command



If a printer is connected to the data interface, pressing the PRINT key (6) initiates the print command.
Note:
In animal weighing mode , pressing this key starts a measuring cycle.

Configuration - file

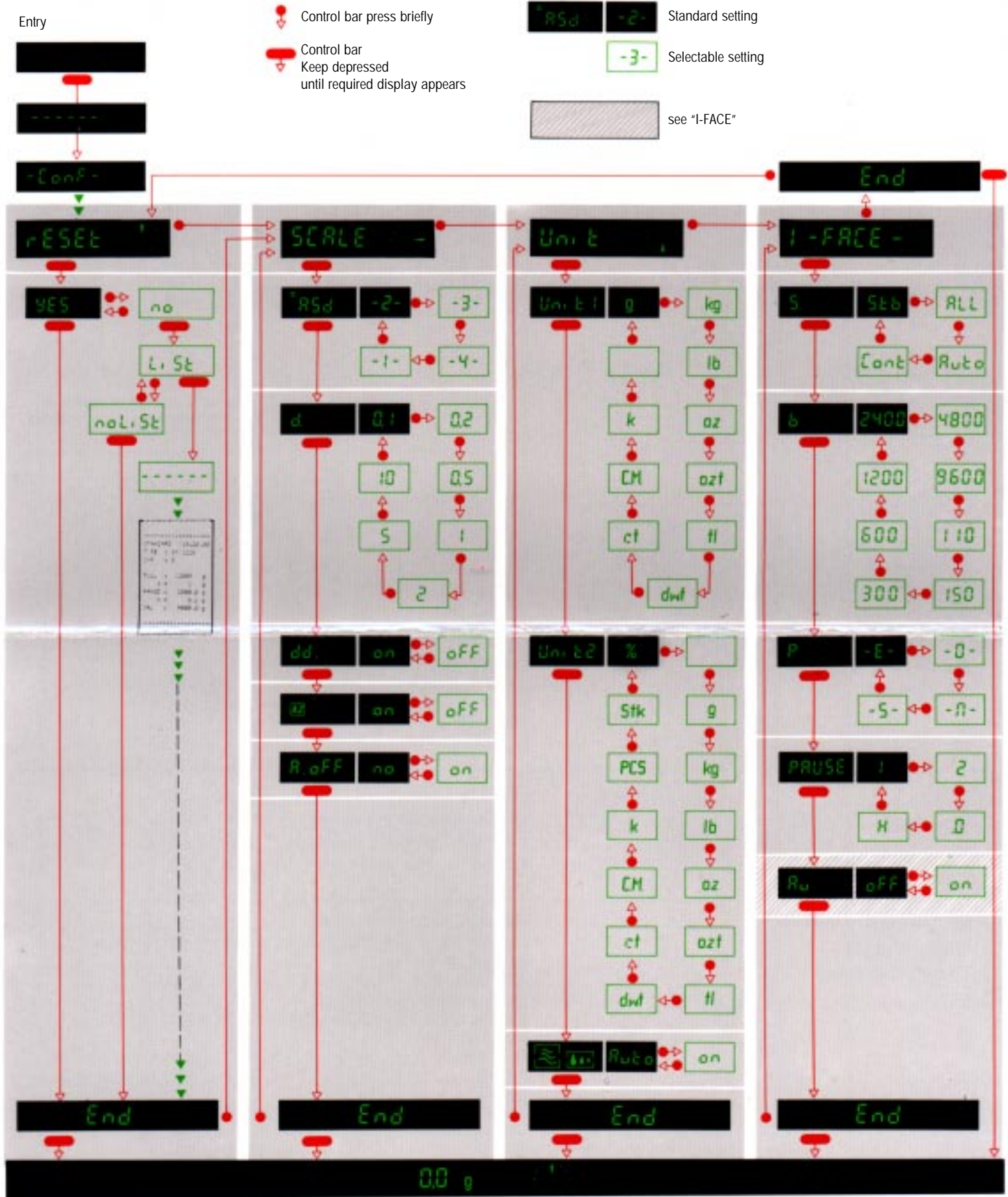

- Important:**
- If you do not press control bar for 40 sec, the scale will automatically return to weighing mode. Modified settings will be recorded.
 - Return to weighing mode, **always** with control bar depressed till zero indication.
 - Display changes automatically.