




## Tour the Controls Palette



The Controls Palette is used to start/stop data collection, set the data monitoring mode, set the sampling rate, delete data runs, define Start/Stop conditions, and playback synchronized video.

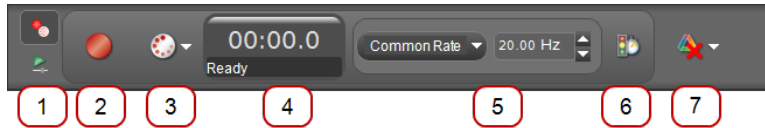





### 1 Record / Play Selection

Icon	Description
Toggle between:	
	Recording Mode Use to switch to recording Controls Palette
	Playback Mode Use to switch to playback Controls Palette

### 2 Record / Monitor

Icon	Description
Capstone toggles between:	
	Record Begin recording all measurements from each connected sensor
	Monitor Start monitoring measurements from connected sensors

### 3 Record Mode Selection

Icon	Description
Toggle between:	
	Continuous Mode Set to record data at the current sample rate settings
	Keep Mode Set to keep data points when commanded
	Fast Monitor Mode Set to acquire data without recording; required for high speed sampling in Scope and FFT displays
 See <a href="#">"How do I record data in PASCO Capstone? ": 78</a>	