

Bio-Plex Pro Mouse Diabetes Assays

Quick Guide

For Use with	Instruction Manual #
Mouse Diabetes Assays	10000092626

This guide can be used to prepare and run a full 1 x 96-well assay plate. For more information on a given step, refer to the complete instruction manual. New users can go to bio-rad.com/bio-plex and download the manual, which includes detailed instructions and a list of kit components.

IMPORTANT! Pay close attention to **vortexing**, **shaking**, and **incubation** instructions. Deviation from the protocol may result in low assay signal and assay variability.










Initial Preparation

1. Plan the plate layout.
2. Start up/warm up the Bio-Plex System (**30 min**).
 - Bring assay buffer, wash buffer, and sample diluent to room temperature (RT). Keep other items on ice until needed
 - Begin to thaw frozen samples
 - Prepare 1x wash buffer. Mix 10x stock by inversion to ensure all salts are in solution. Then dilute 1 part 10x wash buffer (60 ml) with 9 parts distilled water (540 ml)
3. Prime the wash station for flat bottom plate.
4. Calibrate the Bio-Plex System by following the prompts in Bio-Plex Manager Software. This can be done now or during an assay incubation step.
5. Reconstitute a single vial of standards in **500 μ l** of a diluent similar to the final sample type or matrix as shown below. **Vortex** for **5 sec** and incubate **on ice** for **30 min**.

Bio-Plex Pro Mouse Diabetes Assays Quick Guide

Sample Type	Diluent for Standards	Add Bovine Serum Albumin (BSA)
Serum and plasma	Standard diluent	None
Culture media, with serum	Culture media	None
Culture media, serum-free	Culture media	To 0.5% final (w/v)

6. Prepare a fourfold standard dilution series and blank as shown below. **Vortex** for **5 sec** between liquid transfers.

Transfer Volume, μ l	128	50	50	50	50	50	50	50	50	
Reconstituted Standard										
Standard Diluent, μ l	72	150	150	150	150	150	150	150	150	150

7. After thawing samples, prepare as shown below.

Sample Type	Diluent	Add BSA	Sample Dilution
Serum and plasma	Sample diluent	None	Fourfold (1:4)
Culture media, with serum	Culture media	None	Neat to 1:10
Culture media, serum-free	Culture media	To 0.5% final (w/v)	Neat to 1:10

For Adiponectin Assay

Serum and plasma	Serum-based diluent	None	Mouse (1:1,600)
------------------	---------------------	------	-----------------

Note: User-defined validation is required for the use of other dilution factors.

8. **Vortex** the 20x coupled beads for 30 sec and dilute to 1x in Bio-Plex Assay Buffer as shown below. Protect from light.

Number of Wells	20x Beads, μ l	Assay Buffer, μ l	Total Volume, μ l
96	288	5,472	5,760

Running the Assay

Note: Make sure all assay components are at RT before proceeding.

1. **Vortex** the diluted (1x) beads. Add **50 µl** to each well of the assay plate.
2. Wash two times with **100 µl** Bio-Plex Wash Buffer.
3. **Vortex** samples, standards, and blank. Add **50 µl** to each well.
4. Cover plate with sealing tape and protect from light with aluminum foil. Incubate on shaker at **850 ± 50 rpm** for **1 hr** at RT.
5. With 10 min left in the incubation, **vortex** the 20x detection antibodies for 5 sec and quick-spin to collect liquid. Dilute to 1x in detection antibody diluent as shown below.

Number of Wells	20x Detection Ab, µl	HP Detection Ab Diluent, µl	Total Volume, µl
96	150	2,850	3,000

6. Wash the plate three times with **100 µl** wash buffer.
7. **Vortex** the diluted (1x) detection antibodies. Add **25 µl** to each well.
8. Cover and incubate at **850 ± 50 rpm**, as in step 4, for **30 min** at RT. Meanwhile, prepare Bio-Plex Manager Software protocol; enter standard S1 values provided in the assay kit.
9. With 10 min left in the incubation, **vortex** the 100x streptavidin-phycoerythrin (SA-PE) for **5 sec** and quick-spin to collect liquid. Dilute to 1x as shown below and protect from light.

Number of Wells	100x SA-PE, µl	Assay Buffer, µl	Total Volume, µl
96	60	5,940	6,000

10. Wash the plate three times with **100 µl** wash buffer.
11. **Vortex** the diluted (1x) SA-PE. Add **50 µl** to each well.
12. Cover and incubate at **850 ± 50 rpm**, as in step 4, for **10 min** at RT.

Bio-Plex Pro Mouse Diabetes Assays Quick Guide

13. Wash the plate three times with **100 µl** wash buffer.
14. Resuspend beads in **125 µl** assay buffer. Cover plate as in step 4 and shake at **850 ± 50 rpm** for **30 sec**.
15. Remove the sealing tape and **read the plate** using the settings below.

Instrument	RP1 (PMT)	DD Gates	Bead Events
Bio-Plex 200*	High	5,000 (low), 25,000 (high)	50
Bio-Plex 3D*	Enhanced	Select MagPlex Beads	50
Bio-Plex MAGPIX*	N/A, use default instrument settings	N/A, use default instrument settings	Default

* Or similar Luminex system.

Bio-Plex and Bio-Rad are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions.

The Bio-Plex Suspension Array System includes fluorescently labeled microspheres and instrumentation licensed to Bio-Rad Laboratories, Inc. by the Luminex Corporation.

Luminex is a trademark of Luminex Corporation.

All trademarks used herein are the property of their respective owner.

BIO-RAD

**Bio-Rad
Laboratories, Inc.**

*Life Science
Group*

Web site bio-rad.com **USA** 1 800 424 6723 **Australia** 61 2 9914 2800
Austria 43 01 877 89019 **Belgium** 32 03 710 53 00 **Brazil** 55 11 3065 7550
Canada 1 905 364 3435 **China** 86 21 6169 8500
Czech Republic 36 01 459 6192 **Denmark** 45 04 452 10 00
Finland 35 08 980 422 00 **France** 33 01 479 593 00
Germany 49 089 3188 4393 **Hong Kong** 852 2789 3300
Hungary 36 01 459 6190 **India** 91 124 4029300 **Israel** 972 03 963 6050
Italy 39 02 49486600 **Japan** 81 3 6361 7000 **Korea** 82 2 3473 4460
Mexico 52 555 488 7670 **The Netherlands** 310 318 540 666
New Zealand 64 9 415 2280 **Norway** 470 233 841 30 **Poland** 36 01 459 6191
Portugal 351 21 4727717 **Russia** 7 495 721 14 04 **Singapore** 65 6415 3188
South Africa 36 01 459 6193 **Spain** 34 091 49 06 580
Sweden 46 08 555 127 00 **Switzerland** 41 0617 17 9555
Taiwan 886 2 2578 7189 **Thailand** 66 2 651 8311
United Arab Emirates 971 4 8187300 **United Kingdom** 44 01923 47 1301

