# Gel Doc™ EZ Imaging System with Image Lab™ Software

## Version 5.1





#### **Notice**

No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopy, recording, or any information storage or retrieval system, without permission in writing from Bio-Rad.

Bio-Rad reserves the right to modify its products and services at any time. This user guide is subject to change without notice. Although prepared to ensure accuracy, Bio-Rad assumes no liability for errors or omissions, or for any damage resulting from the application or use of this information.

Precision Plus Protein standards are sold under license from Life Technologies Corporation, Carlsbad, CA, for use only by the buyer of the product. The buyer is not authorized to sell or resell this product or its components.

Coomassie is a trademark of BASF Aktiengesellschaft. Coomassie Fluor is a trademark of Invitrogen Corp. Excel, PowerPoint, and Windows are trademarks of Microsoft Corporation. GelGreen and GelRed are trademarks of Biotium, Inc. Krypton is a trademark of Thermo Fisher Scientific Inc. Mac, Mac OS, and Numbers are trademarks of Apple Inc. SYPRO is a trademark of Invitrogen Corporation. Intel and Intel Core are trademarks of Intel Corporation. Mitsubishi is a trademark of Mitsubishi Companies. PulseNet International is a trademark of the Centers for Disease Control and Prevention.

SYBR is a trademark of Life Technologies, Inc. Bio-Rad Laboratories, Inc. is licensed by Life Technologies, Inc. to sell reagents containing SYBR Green I for use in real-time PCR, for research purposes only.

Criterion Stain Free gels are covered by U.S. Pat. No. 7,569,130. Bio-Rad Laboratories, Inc. is licensed by Invitrogen Corporation to sell SYPRO products for research use only under U.S. Patent Number 5,616,502.

Copyright © 2014 Bio-Rad Laboratories, Inc. All rights reserved.

Safety and Regulatory Compliance	9
Safety Use Specifications and Compliance	9
Instrument Safety Warnings	
Notice	
Warranty	
Chapter 1 System Overview	13
Gel Doc EZ Imaging System	15
Technical Specifications	
Image Lab Software Capabilities	
Gel Documentation	
Molecular Weight Assessment	18
Quantification	18
Chapter 2 Setting Up the Instrument with Image Lab Soft	tware19
System Requirements	
Image Lab Security Edition	
Computer Specifications	
Installing Image Lab Software	
Connecting the Imager to the Computer	
Installing the Drivers on a Windows XP Computer	
Installing the Drivers on a Windows 7 Computer	
Configuring the Imager	
Initializing an Additional Tray	37
Initializing an Additional Tray	

	Enabling and Disabling Image Lab Secure Mode	. 49 . 52
	Using Groups on a Local Domain	
Ch	napter 3 Image Lab Software Overview	55
	Interface Overview	. 57
	Main Window	. 57
	Main Toolbar	. 58
	Results Data	. 58
	Display Toolbox	. 58
	Start Page	. 59
	Analysis Tool Box	. 59
	Status Bar	. 61
	Menu Commands	. 62
Ch	napter 4 Acquiring Images	69
	Selecting a Sample Tray	. 70
	Setting Up a Default Protocol	
	Acquisition Settings	
	Analyze Image	
	Regression Methods	
	Output Settings	
	Review Protocol Settings	
	Running a Default Protocol	
	Creating a Multichannel Image from Single Images	
	Setting Up a Custom Protocol	
	Selecting a Custom Protocol	
	Running a Custom Protocol	
	Editing Protocols	

Chapter 5 Viewing Images	89
Displaying Gel Images	90
Display Gel Options	90
Zoom Tools	91
Fit in Window	92
Image Transform	92
Image Colors	94
3-D Projection	95
Image Info	96
Displaying Data	96
Analysis Table Options	
Lane and Band Table Measurement Definitions	99
Volume Measurement Definitions	
Lane Profile	
Standard Curve	
Report	105
Chapter 6 Analyzing Images	
Using Auto Analysis	
Analysis Tool Box Tools	
Image Tools	
Correcting a Slanted Gel	
Cropping a Gel Image	
Inverting Data	
Merging Images	
Lane and Bands Tools	
Detecting Lanes	
Using the All Lanes and Single Lane Tools	
Copying Lanes	
Detecting Bands	
Editing the Detected Bands	
Molecular Weight (MW) Analysis Tools	
Quantity Tools	
Annotation Tools	133
Add Annotations	134

Alignment. Text Properties Color Rotate Volume Tools	135 135 135
Chapter 7 Generating Reports	143
Report	143
Report Settings	144
Print Report	147
Print Report to a PDF File	147
Adjust the Printer Settings	148
Chapter 8 Exporting Results	149
Exporting Gel Images	
Exporting Gel Images for Publication	
Exporting Gel Images for Analysis	
Exporting Gel Images to PulseNet International	
Exporting Lane and Band Tables to Excel	
Exporting Volume Tables to File	
Screenshot Tool Export	
Analysis Table Export	
Chapter 9 Image Lab Logs	155
Image Lab Logs.	
Viewing the Instrument LogViewing the System Log	
Viewing the System Log	
Displaying Log Data	
Displaying Data Columns in Logs	
Filtering Data in Logs	
Collapsing or Expanding Data Rows	
Exporting Logs	
Printing Logs	

Chapter 10 Using the Security Edition	167
21 CFR Part 11	167
Standard Mode versus Secure Mode	168
User Names, Groups, and Roles	168
Role Restrictions	170
Starting Image Lab Security Edition	171
Electronic Records	171
Unsecured Documents	172
Secure Documents	172
Modifying Secure Documents	173
Signing Documents	174
Document Logs	175
Viewing the Document Log	176
Annual dia A. Maintanana and On a ifications	477
Appendix A Maintenance and Specifications	
Cleaning the Sample Trays	
UV-B Fluorescent Lamp Replacement	
Error Messages	179
Appendix B Troubleshooting	181
Appendix C Setting Up Users and Groups	183
Setting Up Image Lab Users and Groups	183
User Accounts	183
User Authentication and Group Membership	
Finding the name of your authentication domain	185
Configuring Users and Groups on a Local Computer	186
Configuring Users and Groups on a Network Domain	190
Password Security	194
Password Policy Setting Examples	196
Account Lockout Policy Setting Examples	196
Auditing Windows Event Logs	196
Miscellaneous Security Measures	197

Appendix D Using Bio-Rad's Stain-Free Technology	199
Stain-Free Workflow	200
Electrophoresis with Stain-Free Gels	
Imaging Gels	201
Imaging Blots	202
Appendix E Regression Calculation Methods	203
Appendix F Mitsubishi P93/P95 Thermal Printer	205
Setting up a Thermal Printer on Windows	
Setting up a Thermal Printer on a Mac	206
Glossary	207

# Safety and Regulatory **Compliance**

The Gel Doc™ EZ imaging system is intended for laboratory use only. To help you make informed decisions about safety, we have provided comprehensive operating procedures and safety information in this manual and on labels affixed to the imager. This information will alert you to any potential hazards. It is the user's responsibility to read and understand the safety information and use it for safe operation of the system.

# Safety Use Specifications and Compliance

The Gel Doc EZ imager is designed and certified to meet EN61010, the internationally accepted electrical safety standard, and EN61326 Class A EMC regulations. Certified products are safe to use when operated in accordance with the instruction manual.

This instrument should not be modified or altered in any way. Modification or alteration of this instrument will:

- Void the manufacturer's warranty
- Void the regulatory certifications

Create a potential safety hazard

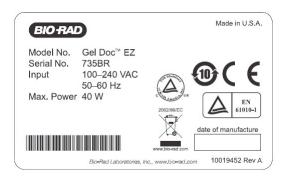


Fig. 1. Gel Doc EZ imager certification label.

For easy customer access, the serial number appears in two locations on your instrument: on the back panel and inside the front door.

# **Instrument Safety Warnings**

Notes, cautions, and warnings are used to highlight certain operating procedures and recommendations. The table below describes how each is used in this document.

Table 1. Notes, Cautions, and Warnings

Icon	Meaning
Note:	<b>Note:</b> A note indicates a special procedure, an exception to normal operation or something else of specific interest to the reader. Notes are preceded by the word <b>Note</b> .
	<b>Caution:</b> A caution precedes an operational step that could damage the instrument or destroy data unless the operator takes certain precautions. Cautions located in the main text are preceded by the word <b>Caution</b> and are accompanied by the Caution symbol in the left margin.

Table 1. Notes, Cautions, and Warnings, continued

#### Icon Meaning

Caution: With the exception of cleaning or replacing light bulbs, refer all servicing to qualified Bio-Rad personnel or their agents. If you experience technical difficulties with the instrument, contact Bio-Rad to schedule service. The instrument should not be modified or altered in any way. Alteration voids the manufacturer's warranty and might create a potential safety hazard for the user.

Caution: If the case interlock is defeated, there is a possibility of UV-B radiation hazard due to UV-B light exposure. Exercise caution when servicing the instrument.

Caution: Bio-Rad is not responsible for any injury or damage caused by the use of this instrument for purposes other than that for which it is intended, or by the modification of this instrument when not performed by qualified Bio-Rad personnel or their agents.

**Caution:** Disconnect the AC power cord before removing the instrument cover.



Warning: A warning precedes an operating procedure that could cause injury to the operator if not followed correctly. Warnings located in the main text are preceded by the word Warning and are accompanied by the Warning symbol in the left margin.

Warning: This instrument must be connected to an appropriate AC voltage outlet that is properly grounded.

## **Notice**

This system is meant for use by specialized personnel who know the health risks associated with reagents normally used in electrophoresis. The UV light source is computer controlled, and proper interlocks are implemented to avoid users' accidental exposure to UV radiation. Bio-Rad Laboratories, Inc. is not responsible for any injury or damage caused by use of this instrument for purposes other than those for which it is intended, or for instrument modifications not performed by Bio-Rad Laboratories, Inc. or an authorized agent.

## **Warranty**

Each Gel Doc EZ imager is protected by a comprehensive instrument warranty agreement. Please read this manual thoroughly so that you fully understand the coverage provided and are aware of your rights and responsibilities. One of the responsibilities of system ownership is regular maintenance. Following the maintenance instructions provided with this manual will help keep your system and peripherals functioning optimally and will protect your investment. Bio-Rad offers a range of comprehensive service agreements that can be tailored to meet your specific needs. Bio-Rad Laboratories is dedicated to your total satisfaction and will be pleased to answer any questions you might have.

# 1 System Overview

This manual describes the Gel Doc™ EZ imaging system, which consists of the Gel Doc EZ imager and Image Lab™ software.



The Gel Doc EZ imager supports multiple applications, including Coomassie and ethidium bromide staining, blue excitation for nondestructive DNA visualization, and stain-free gel imaging.

When coupled with the stain-free gel tray, the Gel Doc EZ imager represents the next generation in stain-free imaging. It features one-button acquisition, yields guick results with higher image quality, and is so simple to use it requires no training, even for users who are not familiar with image analysis systems. The Gel Doc EZ imager's compact size frees up benchspace for running your experiments.

#### 1 | System Overview

The Gel Doc EZ imager is a reproducible and fast label-free SDS-PAGE/native PAGE system that eliminates time-consuming staining and destaining steps. Image Lab image acquisition and analysis software works with the Gel Doc EZ imager to create an automated and time-saving system to image and analyze electrophoresis gels. Data can be viewed, modified, and reported using Image Lab software.

Image Lab protocols can automate gel image acquisition, analysis, and report generation with the push of a single button on the imager.

Researchers can run default protocols or easily design their own custom protocols. Image Lab enables you to view analyzed data, adjust the analysis, and produce customized reports that show precisely the settings that were applied, assuring repeatable results.

# **Gel Doc EZ Imaging System**

The Gel Doc EZ imaging system is a dedicated low-noise detection camera and UV transilluminator-based system with a small footprint. Designed for economy, automation, and ease of use, the user interface consists of a set of four sample trays and a one-touch button to image your gel using a protocol.



Fig. 2. Gel Doc EZ imager, catalog #170-8270.

Note: This catalog number is for the imager with the stain-free tray only. The other trays must be purchased separately.

#### 1 | System Overview

Four sample trays are available for the Gel Doc EZ imager. Each tray supports applications requiring a different set of stains and detection reagents, as shown in Table 2.

**Table 2. Sample Tray Types and Detection Reagents** 

UV Tray Catalog #170-8271	White Light Tray Catalog #170-8272	Blue Tray Catalog #170-8273	Stain-Free Tray Catalog #170-8274
Ethidium bromide	Fast Blast™ DNA stain	SYBR® Green	Stain-free gels
SYBR® Green	Coomassie Blue stain	SYBR® Safe	Stain-free blots
SYBR® Safe	Copper stain	SYBR® Gold	
SYBR® Gold	Zinc stain	GelGreen	
GelGreen	Silver stain		
GelRed			
Flamingo™ fluorescent gel stain			
Oriole™ fluorescent gel stain			
Coomassie Fluor Orange			
SYPRO Ruby			
Krypton			

# **Technical Specifications**

	Hardware Specifications		
Image area	15 x 11.2 cm		
Excitation source	Trans UV B (302 nm peak)		
Detector	Charge-coupled device		
Image pixel size	107.8 x 107.8 μm (in microns)		
Dynamic range	3.0 orders of magnitude		
Pixel density	4,096 gray levels		
Flat fielding	Yes		
Instrument size 43 x 28 x 38 cm			
Instrument weight	Instrument weight 7.3 kg		
Connector descriptions	USB-A connector is used for communication with a PC USB-B connector is used for the instrument interface		
Operating Ranges			
Operating voltage 110/115/230/240 VAC Nominal			
Operating temperature	10–28°C (21°C recommended)		
Operating humidity	< 70% noncondensing		
	Equipment Ratings		
Input voltage range 100–240 VAC			
Input frequency range	50–60 Hz		
Power	40 W		
The Gel Doc EZ imager is	for indoor laboratory use only.		
	Automation Capabilities		
Workflow automated execution	d Controlled by a protocol via setup for gel activation, image capture, analysis, and reporting		
Workflow reproducibility	100% repeatability via recallable protocols, from gel activation and image capture to quantitative analysis and reports		
Autoexposure	2 user-defined modes (intense or faint bands)		

# **Image Lab Software Capabilities**

Image Lab software runs automated protocols on the Gel Doc EZ system for routine gel imaging documentation and analysis.

#### **Gel Documentation**

Gel electrophoresis is a common way to separate, identify, and purify proteins or nucleic acids. The Gel Doc EZ system enables you to image and print gel images for documentation in laboratory notebooks or to export images for publication or presentation.

# **Molecular Weight Assessment**

Sodium dodecyl sulfate polyacrylamide gel electrophoresis (SDS-PAGE) and agarose gel electrophoresis are used to separate proteins or nucleic acids according to their size. Protein or nucleic acid molecular weight standards provide a reference for estimating the molecular weight of the proteins or nucleic acids being tested.

## Quantification

The components in a sample (bands) are quantified to determine relative amounts of each component in a sample or to compare the amount of a sample component relative to a standard. Assessing the purity of a sample requires the quantification of all components of a sample relative to each other. The results are expressed either as a percentage of all bands identified (band%) or as a percentage of all signals in the sample lane (lane%). Determining expression levels of a protein or nucleic acid requires quantification among samples. Data can be reported as either relative values, if the quantity is unknown, or as an absolute value, if a standard of known quantity is present.

# **Setting Up the Instrument** with Image Lab Software

# **System Requirements**

Image Lab™ software runs on Microsoft Windows XP Professional, Microsoft Windows 7, and Mac OS X. Images scanned at high resolution can be quite large. The amount of memory required for using the program is determined by the size of the images you scan and analyze.

For this reason, we recommend that you archive images on a network file server or on removable storage media. Bio-Rad can also provide an appropriate computer to use with this system. Contact your local Bio-Rad representative for more details.

# **Image Lab Security Edition**

Note: The system requirements for Image Lab Security Edition are the same as those for the Standard Edition. The software must be installed on a computer running the Windows XP Professional or Windows 7 operating system to take advantage of the secure mode features.

# **Computer Specifications**

Specifications	Minimum	Recommended
Operating system	Windows XP SP3	Windows XP SP3 Professional
	Windows 7, 32- and 64-bit	Windows 7 Professional, 64-bit
	Mac OS X 10.6	Mac OS X 10.6
Processor	Pentium 4 or equivalent (Windows) at 2.0 GHz	Intel Core 2 Duo 2.0 GHz or higher

#### 2 | Setting Up the Instrument with Image Lab Software

Specifications	Minimum	Recommended
Hard disk space	20 GB	>100 GB
Memory (RAM)	1,024 MB	>1,024 MB
Ports for connecting instrument	1 free USB 2.0 port	1 free USB 2.0 port
Other software (optional)	Microsoft Excel 2003 or later (Windows)	Microsoft Excel 2003 or later (Windows)
	Office 2008 or iWork software (Mac)	Office 2008 or iWork software (Mac)

# **Installing Image Lab Software**

The process of installing Image Lab differs depending on your operating system. This section explains how to install Image Lab software on a Windows PC and on a Mac.

Note: During the Windows installation process, you are prompted to install Image Lab Standard Edition or Security Edition. In order to enable secure mode, you must have Image Lab Administrator role privileges. Bio-Rad recommends that you create the required groups and assign the Image Lab roles to users within those groups before installing Image Lab software.

See User Names, Groups, and Roles on page 168 for additional information about the required Image Lab roles, groups, and users.

See Setting Up Users and Groups on page 183 for information about setting up groups, user names, and passwords.

#### To install Image Lab on a Windows PC

1. Insert the Image Lab software CD in your CD-ROM drive.



The Image Lab installer automatically launches.

- On the Welcome screen, click Next.
- Accept the license agreement and click Next.
- On the Edition Selection screen, choose the edition to install.



Note the following.

- If you are licensed to install the Security Edition but choose to install the Standard Edition, you will need to uninstall the Standard Edition and install the Security Edition before you can use it.
- If you are licensed to install the Security Edition and choose to install it, only a user assigned the Image Lab Administrator role (or group) privileges can enable and disable secure mode.

Note: A user assigned the Image Lab Administrator role will not necessarily also be the network or IT administrator. You must be assigned the correct role within Image Lab to enable or disable secure mode.

For more information, see Setting Up Image Lab Security Edition on page 38.

- If you are not licensed to install the Security Edition but choose to install it, you will be prompted for a license key when you start Image Lab.
- 5. Click Next.

6. A screen appears requesting your license code.

**Note:** Your 18-digit license code can be found in the Image Lab product folder pocket.



Enter the 18-digit code in the three text boxes. The software verifies the code.

**Tip:** If you do not know or do not have access to the code, contact your Bio-Rad customer service representative.


**Note:** Until you provide a license code, Image Lab will function only in standard mode.

- 7. Click Next.
- 8. On the Install Location screen, accept the default location or click Change and browse to another folder.
- 9. Click Next.


10. On the Ready to Install the Program screen, click Install.




The wizard installs Image Lab.

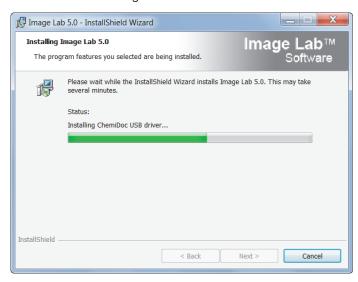

11. When the installation is complete, you are prompted to display the Release Notes and/or the Windows Installer log.

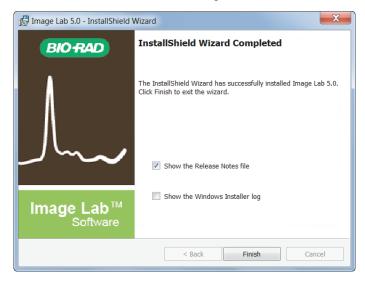

The following screen shot is an example of the Windows Installer log.

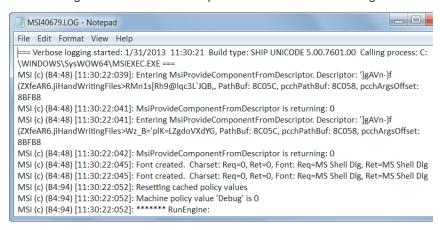

- 12. Select or clear the checkboxes in the Install Wizard Completed dialog.
- 13. Click Finish to exit the wizard.

The Image Lab icon appears on your desktop. Follow the instructions in the next section to connect your system.


#### To install Image Lab on a Mac

- Insert the Image Lab software CD in your CD-ROM drive.
- Double-click the CD icon on your desktop to see the folder contents.
- 3. Double-click the file Image Lab.dmg.
- Drag the Image Lab application icon into the Applications folder.


Follow the instructions in Connecting the Imager to the Computer to connect your system.

# **Connecting the Imager to the Computer**

- Install or upgrade Image Lab before connecting the Gel Doc™ EZ imager to the computer.
- 2. Connect the imager to the Image Lab computer using the USB cable provided with the system.
- Connect the Gel Doc EZ imager to AC power using the power cord provided.

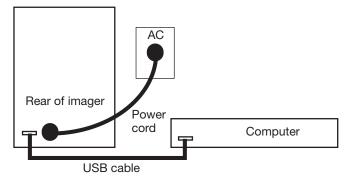

Press the power switch on the side of the imager to turn on power to the instrument.

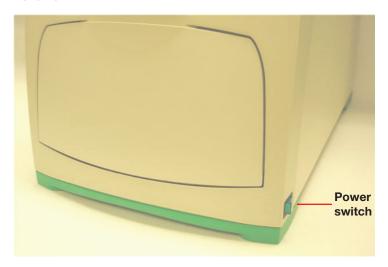

# **Installing the Drivers on a Windows XP Computer**

After the instrument is connected and the power is turned on, the PC automatically recognizes the instrument, and the hardware installation wizard starts, prompting you to install the drivers.


The initial Welcome dialog box appears.



#### Installing the drivers on a Windows XP computer

1. On the Welcome dialog box, click No, not this time, and then click Next.

The dialog box listing the driver that will be installed appears.

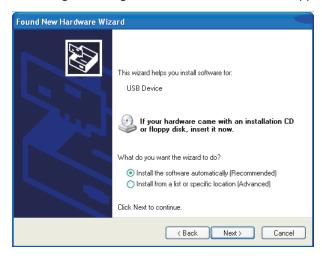

2. By default, the radio button for Install the software automatically (Recommended) is selected.

Keep the default selection and click Next.


The following warning dialog box appears.



#### 3. Click Continue Anyway.

Windows installs the driver.

When the installation is complete, the following dialog box appears.

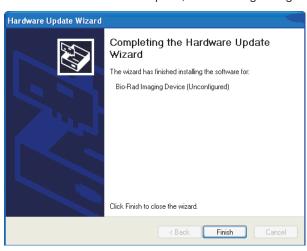

4. Click Finish.

A message appears briefly: Your new hardware is ready to use.


Note: Windows XP might prompt you to go through the driver installation a second time. If this happens, repeat this section until you receive the message "Your new hardware is ready to use."

# **Installing the Drivers on a Windows 7 Computer**

If you are running Windows 7, the device driver is installed during the Image Lab installation process. After successful installation, you see a message similar to the following.



Note: During the installation process you might see a warning similar to the following. You can ignore this warning because it appears even when the driver has been installed correctly.

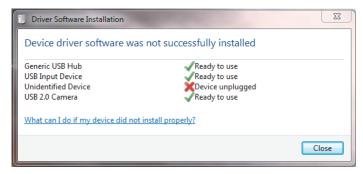

# **Configuring the Imager**

To use the imager with Image Lab software, the imager must first be configured. Imager configuration consists of initializing the sample trays and collecting the dark image.

The imager is configured in the setup wizard, which appears the first time the imager is connected to a computer.

#### To configure the imager

- Turn on the imager and start Image Lab software. The setup wizard appears.
- 2. Complete the steps in the wizard.

Note: You must complete all the steps in the wizard or the imager will not be usable.

The first page of the setup wizard indicates that a new instrument has been found on the system and displays the instrument's serial number.

1. Click Next.




The Dark Image Correction page appears. Correcting the dark image reduces dark current noise generated from the charge-coupled device (CCD). Dark current noise is typical of all CCDs and is a result of the accumulation of charge in the absence of light.

- If a tray is in the imager, remove it and close the tray door.
- Click Acquire Dark Image.

A progress indicator reports the acquisition of the dark image.

When the dark image has been acquired, the wizard prompts you to initialize the UV tray. If you do not have a UV tray, click Skip and go to Step 8.

4. Make sure the UV tray is clean. Then insert the tray into the imager and close the door.

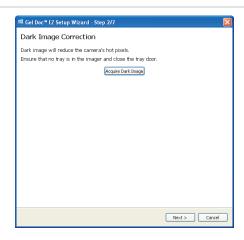



5. Click Next to initialize the UV tray.

A progress indicator reports the initialization of the tray.

When the tray has been initialized, the next screen confirms that the tray has been found and initialized.

- 6. Remove the UV tray from the imager.
- 7. Click Next.

The wizard prompts you to initialize the white tray. If you do not have a white tray, click Next and go to Step 12.

- 8. Make sure the white tray is clean. Then insert the tray into the imager and close the door.
- 9. Click Next to initialize the white tray.


A progress indicator reports the initialization of the tray.





When the tray has been initialized, the next dialog box confirms this and lists the applications you can run with the tray.

- 10. Remove the white tray from the imager.
- 11. Click Next.

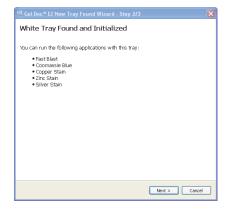

The wizard prompts you to initialize the blue tray. If you do not have a blue tray, click Skip and go to Step 16.

12. Make sure the blue tray is clean. Then insert the tray into the imager and close the door.



13. Click Next to initialize the blue tray.

A progress indicator reports the initialization of the tray.

When the tray has been initialized, the next screen confirms that the tray has been initialized.

- 14. Remove the blue tray from the imager.
- 15. Click Next.

The wizard prompts you to initialize the stain-free tray. If you do not have a stain-free tray, click Skip and go to Step 19.

- 16. Make sure the stain-free tray is clean. Then insert the tray into the imager and close the door.
- 17. Click Next to initialize the stain-free tray.


A progress indicator reports the initialization of the tray.






When the tray has been initialized, the next page confirms this and lists the applications you can run with the tray.

18. Click Next.

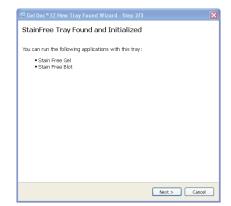

A tray summary page lists all the trays and their status. Trays you initialized appear as ready to use.

19. Click Finish to exit the wizard.

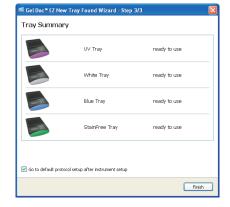

See Chapter 4, Acquiring Images, to configure the default protocols that run when you press the green button on the imager.

# **Initializing an Additional Tray**

After you configure the imager, you can add another tray to the system when you need it.

## To initialize an additional tray

- Insert a new type of tray in the imager. The New Tray Found wizard opens.
- Complete the steps in the wizard.

The wizard identifies the new tray and asks whether you want to initialize it.

1. Click Next.

The wizard begins initializing the tray. A progress indicator reports the initialization of the tray.



When the tray has been initialized, the wizard displays a list of the applications you can run using the tray.

2. Click Next.



The wizard displays the Tray Summary, which shows the highlighted, initialized new trays.

- 3. Do one of the following:
- Click Finish to leave the wizard and display the Default Protocol Setup window, where you can start creating a default protocol for the new tray.
- Clear the checkbox at the bottom of the wizard page to skip setting up a default protocol. Then click Finish to leave the wizard.



# **Setting Up Image Lab Security Edition**

# **Activating and Deactivating Security Edition**

- Before you activate or deactivate Image Lab Security Edition, close any open document files.
- After you change the active status of the Security Edition, you must restart Image Lab.

#### To activate the Security Edition via the Internet

Double-click the Image Lab icon on your desktop to open Image Lab.

The License Code field is populated with the 18-digit license code that you entered when you installed Image Lab.

The Security Edition Activation dialog box appears.

2. Select Activate Via Internet.



Click Activate.



Within about 30 seconds you will receive a confirmation that your Image Lab Security Edition has been activated.



## To activate the Security Edition via email

Double-click the Image Lab icon on your desktop to open Image Lab.

The Security Edition Activation dialog box appears.


The License Code field is populated with the 18-digit license code that you entered when you installed Image Lab.

2. Select Activate Via Create Activation Email.



Click Create Email.

A Save File window appears with the File name field already filled in. Do not change this file name.

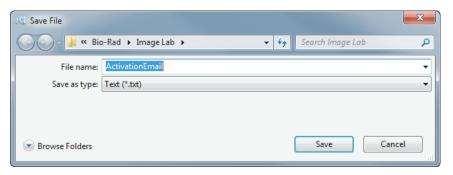

- Click Browse Folders to choose a location for the file, and click Save.
- 5. In your email program, create an email addressed to LSG.TechServ.US@Bio-Rad.com, with the subject line: Request to Activate Image Lab software Security Edition.
- 6. Attach the Activation Email txt file to the email and click Send.

The Bio-Rad Technical Support Department will process your request and reply with an email containing an attachment with your activation code.

- 7. When you receive your reply email, open it and save the attached UnlockCode.txt file to the folder in which you saved the ActivationEmail.txt file.
- 8. On the Security menu, click Activate Security Edition to display the Security Edition Activation dialog box.


9. Select Activate Via Receive Activation Email.




10. Click Receive Email.

An Open File window appears.

11. Navigate to the location where you saved the UnlockCode.txt file, select the file, and click Open at the bottom of the window.

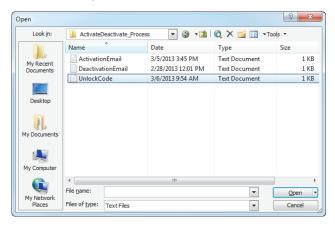

The Security Edition activation complete window appears.

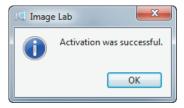

12. Click OK to close the window.

## **Deactivate Image Lab Security Edition**

Image Lab can be installed on more than one computer. Bio-Rad Laboratories recommends that you install the software on only one desktop computer and one laptop. To load the Image Lab Security Edition on a second computer, the software must be deactivated on the first computer before it can be activated on another. You can deactivate the Security Edition automatically via the Internet, or you can deactivate it manually by sending a deactivation email.


### To deactivate the Security Edition automatically via the Internet

- On the Security menu, click Deactivate Security Edition to display the Deactivate Security Edition dialog box.
- 2. Select Deactivate Via Internet.



Click Deactivate. 3.

The system sends a message to the Bio-Rad Technical Support Department, a window stating that your deactivation was successful appears, and Image Lab Security Edition is immediately deactivated.

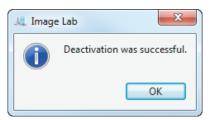


## To deactivate the Security Edition via email

- From the Security menu, click Deactivate Security Edition to display the Deactivate Security Edition dialog box.
- 2. Select Deactivate Via Create Deactivation Email.



Click Deactivate.

A Save File window appears.

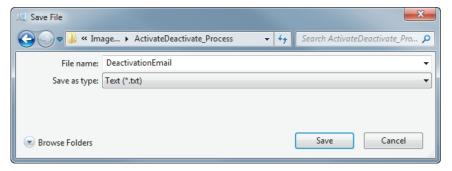

- 4. Navigate to the folder in which you want to save the deactivation email and click Save.
- 5. Create an email addressed to LSG.TechServ.US@Bio-Rad.com, with the subject line: Request to Deactivate Image Lab software Security Edition.
- 6. Attach the DeactivationEmail.txt file to the email and click Send in your email program.


The Bio-Rad Technical Support Department processes your request and deactivates Image Lab Security Edition.

## **Enabling and Disabling Image Lab Secure Mode**


You must have Image Lab Administrator role privileges to switch between Image Lab secure mode and standard mode.

#### To enable secure mode

From the Security menu, click Security Preferences to display the Security Preferences dialog box.

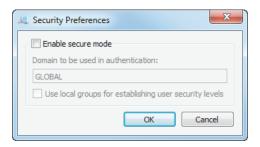

- 2. Select Enable secure mode.
- 3. Click OK to display the Enable Secure Mode authentication dialog box.

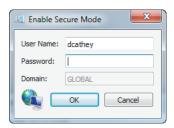


- Enter your Image Lab Administrator user name and password.
- If you are set up on a Windows network server, type the name of the Windows domain in the Domain field.

Note: The default is the domain on which the current Windows user is located, and this name will appear in the field

6. Click OK to save your changes.

A message appears stating that you must restart Image Lab for the new secure mode enable setting to take effect.

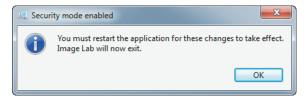


7. Click OK.

The application exits.

Important: For full details on why and how to set your security preferences, see Setting Security Preferences on page 49.

#### To disable secure mode

1. From the Security menu, click Security Preferences to display the Security Preferences dialog box.

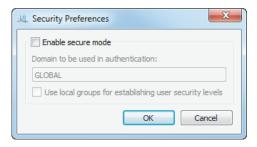

- Clear the Enable secure mode checkbox.
- 3. Click OK.

The Admin Authentication dialog appears.




- Enter your Image Lab Administrator user name and password.
- 5. If you are set up on a Windows network server, type the name of the Windows domain in the Domain field.

**Note:** The default is the domain on which the current Windows user is located, and this name will appear in the field.

6. Click OK to save your changes.



A message appears stating that you must restart Image Lab for the new security settings to take effect.

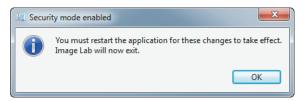

#### 7. Click OK.

The application exits.

# **Setting Security Preferences**

Here are three useful definitions used in these instructions.

- Network domain a remote domain-controlling computer or system which ensures that only authorized users with valid credentials can access and run Image Lab
- Local domain (or local computer) the computer on which Image Lab is running, and which ensures that only authorized users with valid user credentials can access and run Image Lab
- Credentials the valid user name and password that allows or prohibits specific user actions


There are three possible combinations of settings in the Security Preferences dialog box.

### To choose settings in the Security Preferences dialog box

 To set preferences so that only users who are set up on a network domain can use Image Lab: In the Domain to be used in authentication field, enter the name of your network domain. See To find the name of your network domain on page 186 for instructions on how to determine this name.

Note: The default domain name that will appear in this field is the domain on which the current user is logged in.


Do not select the Use local groups for establishing user security levels checkbox.

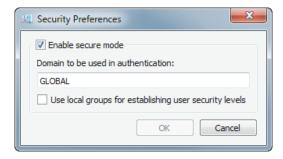

- To set preferences so that only users who are domain users and who are also valid members of specific local groups can run Image Lab:
  - In the Domain to be used in authentication field, enter the name of your network domain. See To find the name of your network domain on page 186 for instructions on how to determine this name.

Note: The default domain name that will appear in this field is the domain on which the current user is logged in.


Select the Use local groups for establishing user security levels checkbox.

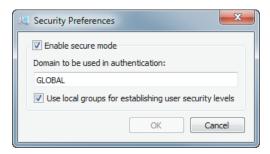

- To set preferences so that only local users can run Image Lab:
  - In the Domain used in authentication field, enter your local computer name. See To find the name of your local domain on page 185 for instructions on how to determine this name.

Note: The default domain name that will appear in this field is the domain on which the current user is logged in.

The Use local groups for establishing user security levels checkbox is grayed out (not accessible).

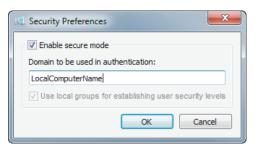

## **Renaming Security Groups**

#### To rename any of the four default Security Groups

From the main menu, select Security > Rename Security Groups.

**Note:** This menu option is visible only if the person logged on to the local computer is logged on as a member of the Windows Administrators group.



- Click in any of the four Group Name fields.
- Enter a new name.
- Click OK to save your changes.

**Note:** The new user group name must comply with standard Windows Local Users and Groups user names rules.

For more information on setting up security groups, see Setting Up Users and Groups on page 183.

# **Using Groups on a Local Domain**

If you choose not to create or use groups on the network domain, set up local groups. Add the authorized users to the groups on the local domain. In the Security Preferences dialog box, select Use local groups for establishing user security levels.

For information about setting up users and groups for Image Lab Security Edition, see Appendix C, Setting Up Users and Groups on page 183.

# **Changing Security Preferences**

Changing the domain that is used to authenticate users is a two-step process. You first authenticate on the first domain, then authenticate on the second domain. This change in domains can be performed in either of two ways. It can be performed by one individual assigned the Image Lab Administrator role on both domains, or it can be performed by two individuals, one with the administrator role on the first domain, and the other with the administrator role on the second domain. See User Authentication and Group Membership on page 183 for more information about using this dialog box.

2 | Setting Up the Instrument with Image Lab Software

# Image Lab Software **Overview**

Image Lab™ image acquisition and analysis software works with the Gel Doc™ EZ imager, stain-free gels, and other gels that are not stain-free to create a reproducible, automated, and time-saving workflow for imaging and analyzing gels.

In Image Lab, a protocol is any combination of imaging, analysis, and report settings that has been saved to run as a single workflow. Researchers can run one protocol repeatedly or easily design a wide range of protocols.

With Image Lab you can view analyzed data, edit the analysis, and produce customized reports that show precisely the settings applied in order to ensure repeatable results.

Image Lab generates two types of files:

- Protocol files describe the parameters for imaging and analyzing your gel images.
- Image files contain the imaged gel, annotations, and analysis performed on the gel. An imaged gel, run according to a protocol file, generates an image file.


Table 3 lists the extensions and icons for the type of files that Image Lab generates.

Table 3. Image Lab file extensions and icons

File type	File Extension	1	Icon	
	Unsigned	Signed	Unsigned	Signed
Protocols	.ptl	.sptl	P	P
Images	.scn	.sscn		

## **Interface Overview**

The following illustration shows the Image Lab main window. This section explains the main software elements.

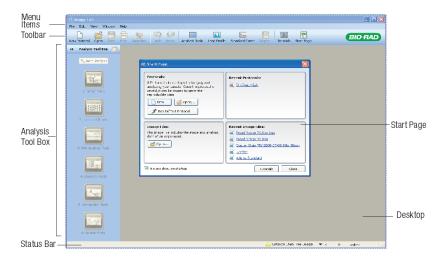

## **Main Window**

Image Lab displays a single main window. All image and protocol dialog boxes that present choices open in the workspace, which is the gray area of the main window.

If many screens are open in the workspace, you can make one active by clicking the title bar at the top of the selected screen. A list of open protocols and image files also appears in the main window menu. Select one to make it active.

You can view complete analyses for images or protocols one at a time or compare image results by arranging screens in the workspace.

## **Main Toolbar**

Many Image Lab tools can be selected by clicking toolbar buttons. The Screenshot tool enables you to send a screen capture of your image to the clipboard or to save it as a file. You can view demonstrations of various functions by clicking Tutorials. The unlimited Undo and Redo buttons enable you to correct missteps easily. The Default Protocols tool at the right end of the toolbar enables you to set up, view, and edit default protocols.




## **Results Data**

Results data associated with gel images can be viewed as an analysis table, a lane profile, a standard curve, or in a report. Different tools for viewing the results data are easily accessible from the main toolbar. These tools are described in Chapter 5, Viewing Images. The views display the analysis for the selected image. All of the views can be displayed at the same time. See Displaying Data on page 96 for details.


# **Display Toolbox**

The display toolbox at the top of every image enables you to display images in the most useful ways. See Chapter 5, Viewing Images for a description of each option.

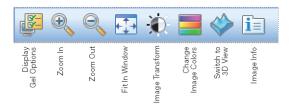


# **Start Page**

The Start Page guides you through creating, opening, and viewing protocols and images.

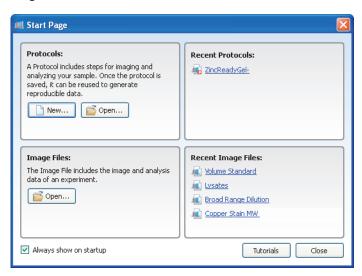

# **Analysis Tool Box**

The Auto-Analysis Auto-Analysis button quickly analyzes images. The remaining tools customize the analyzed data.

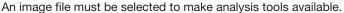

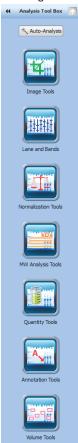


Image Tools enables you to flip, rotate, and crop images and to transform the image files.

Lane and Bands enables you to resize, adjust, and bend lanes and to detect, adjust, add, or delete bands.

Normalization enables you to normalize volume data in multichannel images, so you can correct for sample loading errors in your gels. For more on multichannel images, see the System with Image Lab Software User Guide.

MW Analysis Tools (Molecular Weight Analysis) enables you to choose standard samples, assign standard lanes, and choose a regression method.

Quantity Tools enables you to automatically quantify bands, using either relative or absolute values.

Annotation Tools enables you to add formatted text and arrows to any area of a gel.


Volume Tools enables you to manually quantify an object inside a boundary that you define.

These tools are described in Analysis Tool Box Tools on page 109.

## **Status Bar**

The status bar in the lower right-hand corner of the main window shows the imager and tray in use, and the X and Y values for the cursor position on the image file.

If the default protocol has been set up for the tray in use, its name also appears in the status bar next to the tray name. Finally, the Int status bar displays the intensity value for the image intensity at the cursor position. The maximum data range is 0-4,095, but the range varies depending on the values contained within each image.

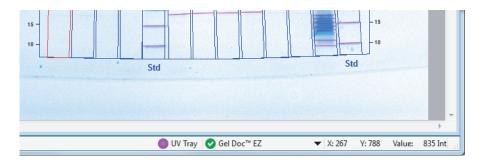

## Menu Commands

The following section describes all menu commands in the File, Edit, View, Window, and Help menus. Many commands are also available on the toolbar or the Start Page.

#### File Menu Commands

**New Protocol** enables you to create a new protocol that contains the necessary steps and choices. See Chapter 4, Acquiring Images for detailed instructions. Protocols can also be altered and stored for reuse.

Open browses the file system to retrieve a previously saved protocol file or image file.

**Recent Images** enables you to open a recent image file.

**Recent Protocols** enables you to open a recent protocol.

**Save** enables you to save a protocol or image file after it is named.

Save As enables you to name and store a protocol or image. Protocols are stored with a .ptl or .sptl extension. Image files are stored with an .scn or .sscn extension.

Close closes the active window.

Close All closes all the windows.

**Export** enables you to export gel images or analysis tables with the following options:

- **Export for Publication** exports a displayed image to a file. You can select from .bmp, .png, .jpg, and .tif formats. The gel displays with any lanes, bands, and annotations that appear on the screen. See Exporting Gel Images for Publication on page 150 for more information.
- Export for Analysis creates a .tif file that retains all gel image data. Analysis data are not included. Use this option to analyze the image in other software such as Quantity One®, FPQuest™, or InfoQuest™FP. See Exporting Gel Images for Publication on page 150 for more information.

- **Export for PulseNet** reduces the image to an 8-bit .tif file. Resolution is limited and file size is restricted to 300 dots per inch (dpi).
- Lane and Band Table to Excel exports your lane and band table data to an Excel (or Numbers on a Mac) spreadsheet.

**Note:** Excel or Numbers must be installed on your computer.

- Lane and Band Table to File exports as a comma-separated values (CSV) file so that your lane and band table can be opened in a database application.
- **Volume Table to Excel** exports your volume table data to an Excel (or Numbers on a Mac) spreadsheet.

**Note:** Excel or Numbers must be installed on your computer.

 Volume Table to File — exports as a CSV file so your volume table can be opened in a database application. See Exporting Volume Tables to File on page 153 for detailed information about exporting files.

See Chapter 8, Exporting Results for more information about exporting files.

Image Info displays information about individual gel and blot images, such as acquisition date and data range, and image capture detail, such as exposure time and illumination source used. Click the Image Details, Analysis Settings, and Notes tabs to display these properties. See Image Info on page 96 for more information.

Page Setup contains print controls such as orientation (landscape or portrait), margins, printer used, and paper size.

Print displays a print preview of the gel and the header information, which includes the filename of the image, the user's name, and the date and time it was printed. The usual Windows Print screen is available as well, enabling you to select a printer and the number of copies to print.

Exit closes Image Lab (after prompting you to save changes to your protocols or images).

#### **Edit Menu Commands**

**Undo** undoes the last action.

**Redo** restores the last action after an Undo.

Screenshot enables you to take a screen shot of the Lane Profile Window, the Standard Curve Window, or the default choice, Current Image View. The screen shot can include the name of the image, and it can be placed on the clipboard or saved in a file.

Default Imager enables users who own two or more imagers to switch between them.

**Instrument Setup** displays information about the instrument, including the instrument name and serial number, camera serial number, firmware version, and the flat field version. The available sample trays are shown with the names of associated default protocols. The dark image time and date are displayed. From this dialog box, you can reset the dark image and the flat field for the white sample tray.

Report Settings enables you to configure reports. This dialog box has three tabs. All of the checkboxes are selected by default. Clear the boxes to exclude information from reports. Your selections apply to all reports until you change them.

- The General tab has options for excluding or reporting information about your gel image.
- The Lane and Band Table tab enables the researcher to choose whether to include all lanes or selected lanes, with appropriate identifiers. Lane profiles can also be included.
- The Volume Table tab enables the researcher to choose appropriate identifiers for the volume table and provides the option of excluding the table from reports.

Preferences enables you to set naming and color preferences for your image files. This dialog box has two tabs.

- The Protocol tab shows presets for naming image files. You can choose to include a designated Prefix, User Name, Date, and/or Time in the name of your image files.
- The Colors tab enables you to choose colors for the graphic elements in your gels, such as Lane Frame, Lane, Band, Band Attribute, and MW Legend. This functionality ensures that these elements are visible. regardless of the color of the gels.

#### View Menu Commands

Image Overview displays the gel image with a red rectangle outlining the area visible in the larger main window. This is useful when you zoom in to a small section of an image.

Image Transform displays a histogram that enables you to adjust the light and dark values of a gel image. This adjustment does not change your data, only the way the data display on your monitor.

**Operations History** displays the sequence of actions performed by both the user and the software.

View System Log displays events related to running Image Lab software, including enabling or disabling secure mode, and the users who log on to or log off of the software.

View (Instrument) Log displays events related to the instrument, including calibrating the instrument and the success or failure of the calibration. This log file is visible only if Image Lab is connected to an instrument.

View (Document) Log displays events related to creating and modifying secure protocol and image files. This log file is created only when Security Edition is enabled.

## **Security Menu Commands**

Note: This menu is only visible in Image Lab Security Edition. For more information about Security Edition, see Chapter 10, Using the Security Edition.

Security Preferences allows the Image Lab Administrator to enable and disable secure mode. In this dialog box, the administrator chooses the domain to be used for authentication and whether to use local groups for security levels.

Rename Security Groups allows a person logged on as a member the Windows Administrators group to change any of the four default Image Lab security group names (TDS\_Administrator, TDS\_User, TDS\_Tech, and TDS\_Guest).

**Note:** This menu option is visible only if the person logged on to the local computer is logged on as a member of the Windows Administrators group. Only users who are logged on as a member of that group maintain the authority to change any of the four default Image Lab security group names. Any changes made to these Security Group names must match the names your Windows system administrator has given those groups.

Sign Document enables users to sign images and protocols. Users enter their user name and password and provide a reason for signing. When the document is signed, the reason is saved in the System Log file.

#### Window Menu Commands

The Window controls enable you to show and hide multiple open image files in your workspace. A list of all currently open images and protocols appears in this menu.

**Tile** aligns all open image files so they are visible at the same time.

**Tile Horizontal** aligns all open image files from top to bottom.

**Tile Vertical** aligns all open image files from left to right.

Cascade stacks all open image files and protocols with overlapping title bars, so each one can be easily chosen for viewing.

Imitate Zoom changes the zoom setting of all open images to the same zoom setting as the current image file.

**Imitate Transform** changes the brightness and contrast of all open images to the same transform settings as the current image file.

Next cycles through all open image files from oldest to newest.

Previous cycles through all open image files from newest to oldest.

## **Help Menu Commands**

Image Lab Help displays the help system.

**User Guide** displays the instruction manual in .pdf form.

**About** displays Image Lab software version and release date.

3 | Image Lab Software Overview

# 4 Acquiring Images

Acquiring an image requires a set of steps in a workflow called a protocol. In Image Lab™ software, a protocol can include any combination of settings for imaging, analyzing, and reporting that has been saved to run as a single workflow.

Users can choose a protocol to be linked to a tray type by default. Such a default protocol is associated with:

- A specific type of tray
- The green Run button on the front of the Gel Doc™ EZ imager
- The user's computer system login ID

The advantage of a default protocol is that it can be run with a single press of the Run button on the imager. Custom protocols are run by clicking a software button in Image Lab software. In all other respects, default protocols and custom protocols are similar.

Typically, you create a default protocol for the applications you run most often with a particular tray. Each user can set up four default protocols (one for each type of sample tray) to run applications. For less frequently run applications, a custom protocol can be created to capture the setup steps for later use.

After it is set up in Image Lab software, a default protocol is run by placing a gel in a sample tray, inserting the sample tray in the imager, and pressing the Run button on the front of the instrument. The imager automatically runs the default protocol assigned to the type of tray inserted. If you have only one tray, it will automatically be considered the default tray.

# **Selecting a Sample Tray**

The following tables list ways of imaging gels stained with various detection agents and which trays to use in each instance. If your stain is not listed here and you are unsure of the appropriate tray, contact Bio-Rad Technical Support for assistance. Stain-free gels and stain-free blots should be imaged in the stain-free tray.

**Table 4. Applications for Nucleic Acid Gels** 

<b>Detection Reagent</b>	Primary Tray	Alternate Tray
Ethidium bromide	UV	None
SYBR® Green	Blue	UV
SYBR® Safe	Blue	UV
SYBR® Gold	Blue	UV
GelGreen	Blue	UV
GelRed	UV	None
Fast Blast™ DNA stain	White	None

**Table 5. Applications for Protein Gels** 

<b>Detection Reagent</b>	Tray
Coomassie Blue stain	White
Copper stain	White
Zinc stain*	White
Flamingo™ fluorescent gel stain	UV
Oriole™ fluorescent gel stain	UV
Silver stain	White
Coomassie Fluor Orange	UV
SYPRO Ruby	UV
Krypton	UV

<sup>\*</sup> Recommended for imaging only.

# **Setting Up a Default Protocol**

Setting up a protocol involves three major processes:

- Acquisition settings
- Analyze image
- Generate output

Each process comprises a set of steps. The appropriate options appear as each protocol step is selected.

Important: You must configure Acquisition Settings (steps 1 through 6 in the following procedure) to run an application. However, it is not necessary to configure Analyze Image and Report Settings unless you want to calculate the molecular weight automatically and generate a report.


## To set up a default protocol

1. Click the Default Protocols button in the toolbar to set up a default protocol.


Caution: If the imager is not connected to your computer, a message appears asking whether you want to run a simulation of the imager. Click No. Turn on the imager. Do not set up a protocol in simulation mode. Protocols set up in simulation mode cannot be run.

#### 4 | Acquiring Images

A Default Protocol Setup window displays each sample tray type as well as buttons that enable you to create a protocol or browse to an existing protocol. Trays that have not been initialized are unavailable for editing or browsing. Note that after you create a default protocol, the Create button changes to View/Edit so you can review or change the protocol settings.

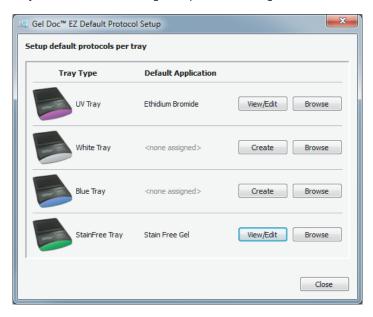

2. Click Create for the sample tray type for which you want to create a protocol. The first Protocol Setup window appears. The title bar identifies the protocol as a default protocol. The imager type and sample tray type are also shown.

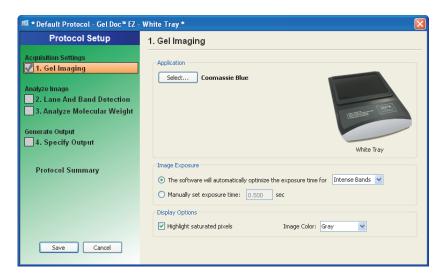

The three major processes are listed in the left pane. Numbered major steps in each process appear under these headings. To select options under a protocol step, select the accompanying checkbox. Options appear in the right side of the window. To disable any step, clear its checkbox.

## **Acquisition Settings**


## Step 1. Gel Imaging

The Gel Imaging checkbox is already selected. The right side of the Protocol Setup window displays Gel Imaging options.

## **Application Options**

### 1. Select an application

Click Select and choose an application from the dropdown menus.

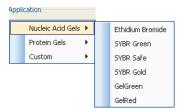

The menus list only applications appropriate for the selected sample tray. When you choose an application, any required filter or illumination source displays in the Protocol Setup window.

**Note:** Custom applications are available only for custom protocols. See Setting Up a Custom Protocol on page 84 for more information about custom applications.

#### 2. Set the image exposure

If you select the first image exposure option, Image Lab acquires an image using the optimal exposure time. This option ensures the best use of the dynamic range. You can automatically optimize for intense or faint bands.



If you choose Intense Bands from the menu, exposure is optimized for all bands.

- If you choose Faint Bands, exposure time is longer, making faint bands more visible. More prominent bands might be overexposed.
- Select manual exposure to override automated imaging. Exposure time can range from 0.001 to 10 sec.

After a gel is imaged with automatic exposure optimization, the previous exposure time displays in the protocol as a guide so you can manually adjust the exposure time if necessary.



You can also view the exposure time of the image later, in the Image Info window (see Image Info on page 96 for more information).

### Set the Display Options

**Highlight saturated pixels** — select the Highlight saturated pixels checkbox to see any saturated pixels in red, which shows how much of the gel image is saturated. Saturated pixels are beyond the maximum quantifiable range of the imaging system. This option can be changed later by selecting View > Image Transform.


**Image Color** — select color choice to display the sample image. Viewing the image with a different color scheme can make it easier to see all of the elements. See Image Colors on page 94 for the color choices.

## **Analyze Image**

## Step 2. Detect Lanes and Bands

To analyze the gel automatically, Image Lab must detect lanes and bands on the image. Lanes are detected automatically, and then the background is automatically subtracted. Refer to Lane and Bands Tools on page 113 for details. You can customize band detection with the following options.

Select the Lane and Band Detection checkbox in the left pane of the Protocol Setup window.

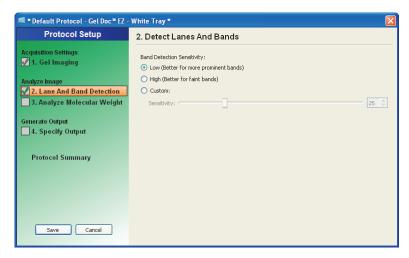

The Protocol Setup window displays lane and band detection options:


- **Low Band Detection Sensitivity** this option sets detection at a low level for images with prominent bands. Faint bands are not detected.
- **High Band Detection Sensitivity** this option sets detection at a higher level for images that are faint. Extraneous bands can be removed later using the Band Tools in the Analysis Tool Box. See Lane and Bands Tools on page 113.

**Custom** — select a numerical value between 1 and 100 to choose the best detection sensitivity for your sample.

Note: When low or high band detection sensitivity is used, these numerical values are set: low sensitivity = 25; high sensitivity = 75.

## Step 3. Analyze Molecular Weight

- 1. To analyze molecular weight automatically, select the Analyze Molecular Weight checkbox in the left pane of the Protocol Setup window. When this checkbox is selected, the software calculates the molecular weight for each band based on the specified standard.
- 2. To estimate the size of the molecules in the bands of your gel, enter the standards you are using and specify the lanes in which the standards are placed.

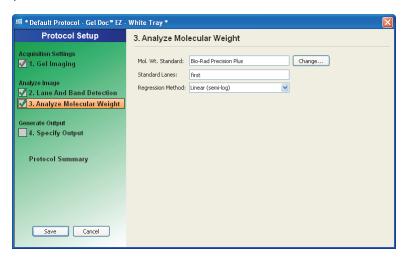


### **Standards**

Determining molecular weight depends upon selection of the proper protein standards. Many protein standards are available from Bio-Rad, including:

- Precision Plus Protein<sup>™</sup> standard
- Prestained SDS-PAGE standards, broad range
- Prestained SDS-PAGE standards, low range
- Prestained SDS-PAGE standards, high range

Bio-Rad standards cannot be edited. However, these standards can be copied, and the copies can be renamed and edited.

The Change button opens the Manage Standards dialog box, in which you can add third-party standard samples and edit your list.

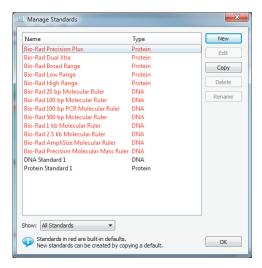

#### Standard Lanes

Choose the lane that contains your standards by typing lane numbers or the words First and Last in the Standard Lanes field. The format is xx, xx, xx, ... where xx is the lane number. For example, if you run an 18-well gel and you want your standards in lanes 1, 10, and 18, enter First, 10, and Last.

Note: Lane detection works best when standards are placed in the first and last lanes.

## **Regression Methods**

A regression method is used to calculate the molecular weight of the unknown bands. The software uses the relative front and molecular weight values of the standard bands to calculate the standard curve. The standard curve is then used to calculate the values of the unknown bands. The shape of the standard curve is based on the selected regression method. Choose one of four regression methods.

The methods and the minimum number of standard bands required for each type are shown in the following table.

Table 6. Regression methods and minimum number of bands

Regression Method	Minimum Number of Standard Bands
Linear (semilog)	2
Point-to-point (semilog)	2
Logistic	5
Cubic spline	5

If you do not have enough data points for the selected method, the molecular weight of the unknown bands is not calculated.

Select the appropriate method for the gel type:

Gradient gels: The linear (semilog) regression method works well for these gels because the mobility of the bands is linear to the log of their molecular

#### 4 | Acquiring Images

weight. As an alternative, the point-to-point (semilog) method can be used if the R<sup>2</sup> value is not sufficient.

Fixed percentage gels: These gels have a nonlinear relationship between the mobility and the molecular weight. For these gels, choose the logistic or cubic spline regression method.

You can check how well each regression method fits the data in the standard curve window (for more information, see Standard Curve on page 104). The linear (semilog) regression method provides a measurement that describes how well the standard curve fits the data R<sup>2</sup> value. The closer the R<sup>2</sup> value is to 1.0, the better the data fit the standard curve

The molecular weight of each band is displayed in the Analysis Table in the Mol. Wt. column. Refer to Molecular Weight (MW) Analysis Tools on page 123 for more information about molecular weight.

For information about the calculations behind the regression methods, see Appendix E, Regression Calculation Methods.

## **Output Settings**

Image Lab prints to the default printer unless you select otherwise.

You have three choices for specifying output:


- You can automatically print the image. This is useful for gel documentation, especially if a thermal printer is connected. See Appendix F, Mitsubishi P93/P95 Thermal Printer for instructions on setting up the Mitsubishi P93/ P95D thermal printer
- You can automatically print the report, including all accompanying analysis information.

**Note:** You cannot print a report on a thermal printer.

You can display the image on your computer.

## Step 4. Specify Output

Select the Specify Output checkbox in the left pane of the Protocol Setup window to display the output options.

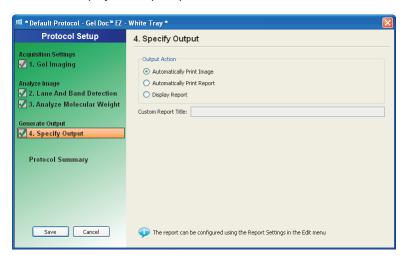

In the right pane, you can choose whether to automatically display or print an image or a report after a protocol runs.


The report can be customized with the options in Edit > Report Settings.

For more information about reporting options, see Chapter 7, Generating Reports.

## **Review Protocol Settings**

### **Protocol Summary**

 Click Protocol Summary in the left pane of any Protocol Setup window to see a review of all protocol settings.

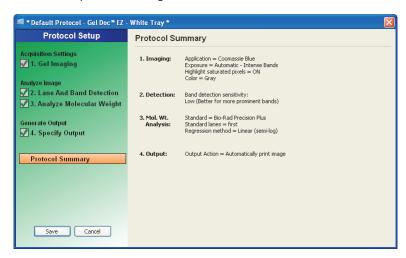


2. Click Save in the left pane to save the default protocol.

## **Running a Default Protocol**

#### To run a default protocol

- 1. Place a gel on the appropriate sample tray and insert the tray into the imager until the magnet grabs the tray.
- 2. Close the door.
- 3. Press the green Run button on the front of the imager. The default protocol runs automatically.

The default protocol appears in Image Lab with a green background. After the protocol runs, the resulting image appears and the protocol window remains open behind it. The application name appears in the status bar below the gel preview window.

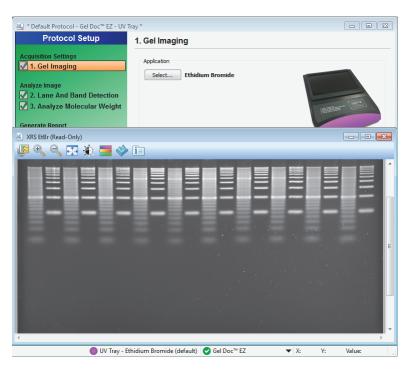

## **Creating a Multichannel Image from Single Images**


You can create a multichannel image from existing single images or from single channels in other multichannel images. Only images with the same aspect ratio can be combined in a multichannel image. These images are not linked to one another. Therefore, when you make a change to one image, it is not propagated to the other images. For information on how to create a multichannel image from single channel images, see the chapter on acquiring images in the ChemiDoc™ MP System with Image Lab Software User Guide.

# **Setting Up a Custom Protocol**

Creating a custom protocol is similar to creating a default protocol.

### To create a custom protocol


- 1. Choose New in the Protocols box on the Start Page or click New Protocol in the toolbar. The Protocol Setup window appears.
- 2. Select an application from the Application dropdown menus.

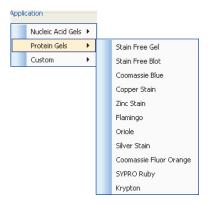

The menus list all predefined applications you can run on the Gel Doc EZ imager.

**Note:** If you want to image a gel using a detection reagent not listed in the menus and do not find an equivalent reagent in the menus, contact Bio-Rad Technical Support for protocol recommendations.

Choose Custom to run an existing application with a new name or when you have an application unlike any existing applications. Selecting Custom > Manage Custom Applications in the Application menu displays the Manage Custom Applications dialog box. Custom applications you save are listed in this dialog box. An Edit button enables you to make changes to custom applications.

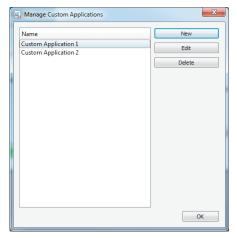


Note: You cannot create custom applications for the stain-free tray.

- To create a custom application, click New. The Create Custom Applications dialog box appears.
- 4. Enter a unique application name.
- 5. The remainder of the setup instructions are the same for both default and custom profiles. Go to step 2, Set the image exposure on page 74 in Setting Up a Default Protocol and follow the instructions.

## **Selecting a Custom Protocol**

Use the same protocol repeatedly by clicking Open in the Protocols box on the Start Page or by clicking Open in the toolbar. You can also select recently used protocols or image files from the lists on the Start Page.

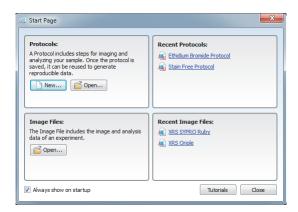

## **Running a Custom Protocol**

#### To run a custom protocol

- Open a protocol or create a new one.
- Place your gel on the appropriate sample tray and insert the tray into the imager until the magnet grabs the tray.
- Close the door.
- Click Run Protocol in the left pane of the Protocol Setup window. The custom
  protocol runs automatically.

After the protocol runs, the resulting image appears and the Protocol window remains open behind it. The application name appears at the bottom of the gel preview window.

# **Editing Protocols**

Protocols can be changed or renamed using the many tools in Image Lab.

## To edit protocols

1. Choose File > Open to open a saved protocol.

Alternatively, you can click Open in the Protocols box on the Start Page. You see the same set of dialog boxes and choices described in the section, Setting Up a Custom Protocol on page 84.

- 2. To replace an existing protocol, edit the protocol and save your changes.
- 3. To create another protocol, save the edited protocol with a different name.


4 | Acquiring Images

# 5 Viewing Images

After a gel is imaged, the image appears in the workspace. Many controls are available to optimize viewing and to analyze the image.

The following screen shot shows a gel image with band and lane detection as well as annotations. The labels are overlays that you can display or hide.

There are many ways to view the data associated with the results. You can view data as an analysis table, a standard curve, and a report.

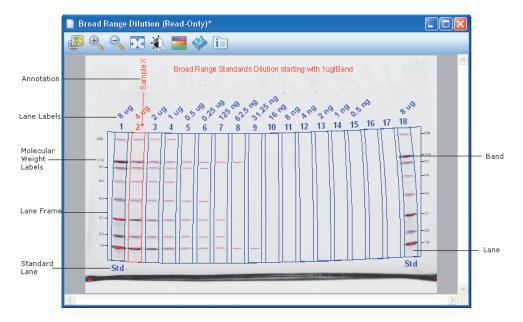

# **Displaying Gel Images**




The display toolbar is located above the gel image. Each tool is described in the following sections.

# Display Gel Options



This section describes the settings in the Display Gel Options dialog box.

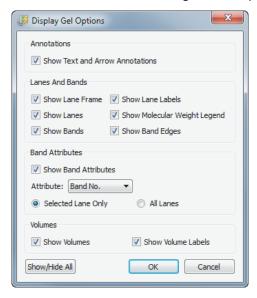

#### **Annotations**

You can choose whether to show text and arrow annotations that have been drawn on the image.

### **Lanes and Bands**

You can turn on or off any image overlays, such as lane frames, lanes, bands, lane labels, molecular weight legends, and band edges.

#### **Band Attributes**

You can show the following attributes for selected lanes or for all lanes.

- Band number
- Band label
- Molecular weight
- Relative front
- Volume
- Absolute Quantity
- Relative Quantity
- Band %
- Lane %

### **Volumes**

If you drew volume boundaries on the gel, you can display the boundaries and their volume labels.

# Zoom Tools



The zoom tools resize the gel image. Click the magnifying glass with the plus sign to make the image larger. Click the magnifying glass with the minus sign to make the image smaller.

You can also zoom in on an area using the right mouse button. Right-click and drag to select the area you want to magnify. You can also resize the image by rightclicking and using the scroll wheel on your mouse.

Tip: You can return to the original view by right-clicking anywhere on the image.


# Fit in Window

If you zoomed in on an area of an image, this button brings the entire image back into view.

# **Image Transform**



The Image Transform dialog box adjusts image brightness and contrast, optimizing the image display so faint details can be seen.

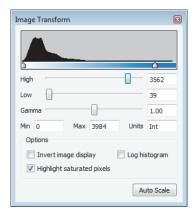

The minimum to maximum range varies depending on the light and dark values present in the image. These adjustments do not change the data. They change only the way the data are displayed. The human eye cannot see as great a range as the image contains.

The frequency distribution histogram shows the total data range in the image and the amount of data at each point in the range.

Auto Scale determines an optimal setting for the image automatically. The lightest part of the image is set to the minimum intensity, and the darkest is set to the maximum.

The High progress indicator determines which intensity value is shown at the maximum gray scale (or other color) in the gel image.

- The Low progress indicator determines which intensity value is shown at the minimum gray scale (or other color) in the gel image.
- The Gamma progress indicator changes the gray scale curve. A value of 1 is linear. A value <1 redistributes a greater proportion of the gray scale to the first half of the intensity values. A value >1 redistributes a greater proportion of the gray scale to the second half of the intensity values.

You can also type numerical values in the boxes next to the progress indicators. Clicking anywhere on the progress indicator bars moves the progress indicator incrementally.

### **Options:**

- Invert image display inverts dark bands on a light background to light bands on a dark background. Light bands on a dark background are inverted to dark bands on a light background.
- **Highlight saturated pixels** when this checkbox is selected, areas of the image with saturated signal intensity (higher than a measurable range) are highlighted in red.
- **Linear or logarithmic histogram** this adjustment changes the y-axis on the histogram to display the number of pixels at each intensity value using either a linear or a logarithmic scale.

# Image Colors **=**




You can choose a color map for your image results file. Viewing the image with a different color scheme can make it easier to see all of the elements in the image, but it does not change your data.

The first eight color choices imitate the colors of stained gels. The remaining choices supply enough color variation to highlight small differences in the image data. The available colors include:

- Gray
- EtBr (ethidium bromide)
- Coomassie
- Stain-free
- SYBR® Green
- SYPRO Ruby
- Flamingo
- Silver

- False color
- Spectrum
- Gold-Silver
- Pseudo

# 3-D Projection

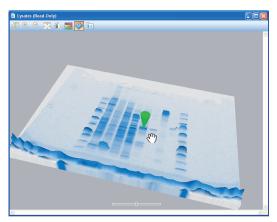

The 3-D view transforms the gel image into a solid three-dimensional model spinning in space with x, y, and z dimensions. Accentuate or diminish the relative heights of data points by pulling the slider at the bottom of the window to the right or left.

### To view the intensity of various bands

- 1. Select the 3-D button in the display toolbar.
- Click and drag the model to rotate it into your preferred view.
- 3. Bring the window into focus by clicking the image.
- 4. Press "C" to display an inverted green cone, which can be dragged around to evaluate the intensity of various bands.
- 5. Press "C" again to hide the tool.

# Image Info

The Image Info dialog box provides information about the active image.

The dialog box has three tabs.

- **Image Details** acquisition and image information appear in this tab.
- **Analysis Settings** settings that were used when the gel was analyzed are displayed here. For example, Band Detection and Molecular Weight Analysis will appear if they were performed.
- **Notes** on this tab you can add notes, point out the types of samples used, and add any other information about the results. You can create custom labels for the lanes in your image.


## **Displaying Data**

You can view results from analyzed data associated with a gel image using an analysis table, a lane profile, a standard curve, or a report.

The buttons in the main toolbar turn these views on or off. You can view your data with all views simultaneously.

# Analysis Table Options

Numerical data associated with an analysis can be viewed in an analysis table. Data from the Lane and Band analysis can be viewed in the Lane and Band Table tab. If volume analysis was performed, these data can be viewed in the Volume Table tab. The buttons above the table provide options for displaying and exporting analysis table data.

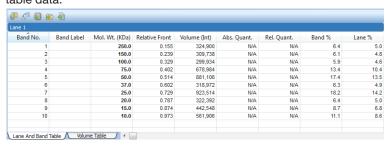

### To change the size of the Analysis Table window

Move your cursor to the top of the window until it changes into a double-headed arrow. Click and drag the edge of the window until you can see all of the data.

**Note:** Resizing the Analysis Table window is restricted when a protocol window is open.

# Display Data Options

The Display Data Options dialog box has three tabs.

Measurements — enables you to choose the measurements to display in the table. Use the arrows to move the columns between the Not Displayed pane and the Displayed pane. By default, all measurements are displayed in the Analysis table.

**Note:** For a description of each Lane and Band measurement type, see Lane and Band Table Measurement Definitions on page 99. For a description of each Volume measurement type, see Volume Measurement

#### Definitions on page 100.

- **Display** enables you to set the display for the analysis table. The following settings appear on the Display tab:
  - **Default display settings** The Move selected lane to top checkbox is selected by default. When you click a lane on the image, the Analysis table scrolls so that the data for the selected lane appears first in either the vertical or horizontal view of the table.
  - □ Per Measurement Precision set the precision (decimal places) for the measurements in the Lane and Band table and the Volume table.
  - □ **Example** shows an example of how measurements will display with the selected measurement and precision settings.
- **Export** enables you to choose how to export the analysis data. The following settings appear on the Export tab:
  - **Export formatting** select checkboxes to include lane headers (Lane and Band table tab only) and/or column headers in the exported file.
  - **Export delimiter** select a delimiter option for the exported file.
    - Comma delimited
    - Tab delimited
    - Use other delimiter (user defined)

## Change Analysis Table Orientation



This button toggles between two table orientations.

Horizontal — displays the lanes/volumes beside each other, so you can scroll through the table from left to right.

**Vertical** — displays the lanes/volumes on top of each other, so you can scroll through the table from top to bottom.

## Copy Analysis Table to the Clipboard

Copies the table data to the clipboard so that you can paste the data into word processing or presentation applications.

**Tip:** It is best to use the vertical table orientation when copying to an 8.5 x 11-inch page, to give the columns enough room to display.

### Export Analysis Table to a File



Exports the table data as a CSV file so you can open it in a database application.

### Export Analysis Table to Excel



Exports the table data to a spreadsheet so you can use the sorting and formula functions to manipulate your data. If you have Excel (PC or Mac) or Numbers (Mac) installed on your computer, the program opens with your spreadsheet displayed.

## Lane and Band Table Measurement Definitions

This section defines the measurements that display in the Lane and Band tab in the Analysis table. Use the Data display options button to choose the columns to display.

**Band Number** — each band in a lane has a unique number, sorted from top to bottom.

Band Label — you can assign a custom label to each band by clicking the Band Label field of the Lane and Band table.

**Molecular Weight** — the molecular weight of the band is calculated based on the user-defined standard and regression method. Italic values indicate extrapolated values. When using nucleic acid gels, the size of the band is displayed in base pairs.

**Relative Front** — values between 0–1 indicate the relative movement of the band from top to bottom.

**Volume** — the sum of all intensities within the band boundaries.

**Abs. Quant.** — absolute quantification of the band.

**Rel. Quant.** — relative quantification of the band compared to the reference band.

**Band** % — percentage of the band's volume compared to all band volumes in the lane.

Lane % — percentage of the band's volume compared to the entire volume of the lane.

### **Volume Measurement Definitions**

This section defines the measurements that display in the Volume tab in the Analysis table. Use the Display Data Options button to choose the columns to display.

**Volume Number** — a unique number is assigned to each volume.

**Volume Label** — software-generated labels for different types of volumes (U – unknown, B – background, S – standard). Label can be changed in Volume Properties.

**Volume** — the sum of all the intensities within the band boundaries.

**Adjusted Volume** — the background-adjusted volume.

**Mean Background** — the mean value of the background.

**Absolute Quantity Volume** — the quantity of the volume based on the standard volumes and the regression method.

**Relative Quantity Volume** — the ratio of the adjusted volume and the adjusted volume of the reference volume.

# **Pixels** — number of pixels inside the volume boundary.

**Minimum Value** — intensity of the pixel with the minimum intensity inside the volume.


**Maximum Value** — intensity of the pixel with the maximum intensity inside the volume.

**Mean Value** — mean value of all pixels inside the volume boundary.

**Standard Deviation** — standard deviation of all pixels inside the volume boundary.

**Area** — area of the volume in mm<sup>2</sup>.



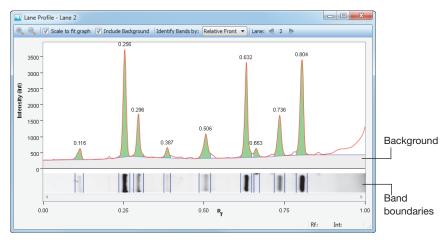

The Lane Profile option shows a cross-section view of a single lane rotated 90°.

### To explore the Lane Profile window

- Use the Next and Previous buttons at the top of the window to page through profiles of the lanes in your image.
- Click on a lane in your image to see its profile.

The title bar identifies which lane profile is in view (Lane 1, Lane 2, and so on).

There are several settings in the title bar.

- Scale to fit graph
- Include Background
- Identify Bands by

These settings, as well as the zoom tools, are global. They apply to all the profiles.

The zoom tools work the same way they work elsewhere in the software. For more information, see Zoom Tools on page 91.

In addition to the graph of the lane intensities, the Lane Profile tool also shows an image of the selected lane below the graph. The transform and color map are applied to the gray-scale image.

As you move your cursor over the profile, the current relative front (Rf) value and the average intensity (Int) value at the Rf value display in the lower-right corner of the Lane Profile window.

### Scale to Fit Graph

You can choose the highest point of the display to define the range of the graph. This provides the best view of the lane profile.

You can clear the Scale to Fit Graph checkbox to display the entire range of possible intensity values in the graph. Doing so allows valid comparisons between different lanes.

### Include Background

When the Include Background checkbox is selected, the Lane Profile window shows the subtracted background under the blue line. The area used for band quantification appears in green under the red line.

When you clear the Include Background checkbox, the area of the lane profile that represents the background of the image does not display.

## Identify Bands by

You can change how the bands are labeled by choosing from the options in the Identify Bands by dropdown list. By default, the bands are labeled by band number.

You can display one of the following attributes:

- Band Number (Band No.)
- Band Label
- Molecular Weight (Mol. Wt.)

- Relative Front
- Volume
- Absolute Quantity (Abs. Quant.)
- Relative Quantity (Rel. Quant.)
- Band %
- Lane %

## **Adjusting Band Boundaries**


Below the profile of each lane there is a strip that displays the bands. Each band is surrounded by a pair of vertical lines delimiting its boundaries. You can move the vertical lines and change these boundaries.

## To change the boundaries of a band

- 1. Hold your cursor over one of the boundary lines until a double arrow appears.
- 2. Click and drag the cursor to the new position.

Note: Boundary lines cannot overlap one another. Therefore, you will not be able to move any boundary line beyond the next boundary line.

# **Standard Curve**

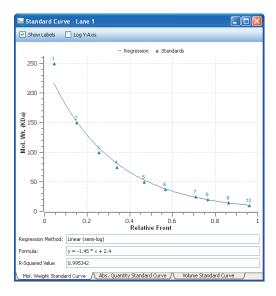

The Standard Curve dialog box displays the best curve fit for the defined standards and the bands relative to this curve for the lane selected in the image. The tabs at the bottom of the dialog box display the standard curves for three different analyses.

Standards appear in green. Unknown bands appear in red. You can toggle the molecular weight display on the y-axis between linear and log scale by clicking the Log y-axis box at the upper left. The regression method you chose in Molecular Weight Analysis Tools appears, as well as the formula (if applicable) and the R<sup>2</sup> value of the regression method.

Tabs in this window enable you to view the molecular weight standard curve, the absolute quantity standard curve, or the volume standard curve.



See Generating Reports on page 143, for information about reports.

## **Analyzing Images** 6

Analysis Tool Box tools are enabled when an image file is opened and in focus. An active or "in focus" window has a darker blue menu bar on a Windows PC. On a Macintosh, the window control icons display more brightly when a window is active. This distinction helps you to identify the active window among many open image files in your workspace.


## **Using Auto Analysis**

Auto analysis detects the lanes and bands and, optionally, calculates the molecular weight of the bands in your image. Click Auto-Analysis in the Analysis Tool Box to do the following:

- Analyze images obtained with protocols that did not include steps for auto detection and analysis
- Change your analysis parameters to reanalyze your images

Note: If you change any settings for an analyzed gel, the initial analysis is overwritten. To preserve both analyses, save each image file with a different name.

### **Auto Detection Settings**

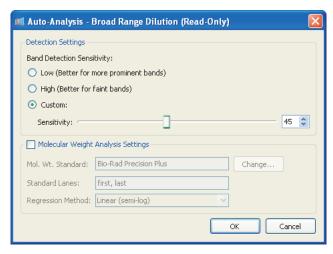

The band detection sensitivity options are as follows:

**Low Band Detection Sensitivity** — sets detection at a low level for images with prominent bands. Faint bands are not detected with this setting.

**High Band Detection Sensitivity** — sets detection at a higher level for images that are faint. Extraneous bands can be removed using the Band Tools in the Analysis Tool Box. See Lane and Bands Tools on page 113.

**Custom** — enables you to set a value between 1 and 100 to select the best detection sensitivity for your sample. You can also drag the sliding bar left or right to set the value.

When Low Band Detection Sensitivity is used, the numerical value is set at 25; when High Band Detection Sensitivity is used, the value is set at 75.

## **Molecular Weight Analysis Settings**

**Molecular Weight Standard** — choose any of the many Bio-Rad standards or another standard you added to your standards list.

**Standard Lanes** — choose or change the lanes in which the standards are placed.

**Regression Method** — four regression methods are available. For more information, see Regression Calculation Methods on page 203.

# **Analysis Tool Box Tools**

All Analysis Tool Box tools customize the analyzed data in image files. These tools are available only when an image file is open. Click a specific image to select among the many windows that might be open in your workspace.

Note: Some tools delete the existing analysis.

#### To access a tool

Click any of the toolbox buttons.

#### To return to the Analysis Tool Box menu

Click the green Up arrow to the right of the tool name.






The image tools enable you to manipulate your images.

#### To display the image tools menu

Click Image Tools.

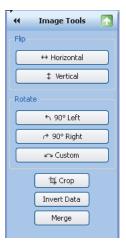

- **Flip** you can flip the gel image horizontally or vertically.
- **Rotate** you can rotate the gel image 90° using the Left or Right buttons. Or you can set a custom rotation using the Custom button.
- **Crop** you can trim the outer edges of your image to any shape or area.
- **Invert Data** you can toggle the image data from positive to negative.
- Merge you can merge a chemiluminescent blot image with a colorimetric image of the same blot.

The sections that follow describe how to use these tools in greater detail.

# **Correcting a Slanted Gel**

#### To correct a slanted gel

- In the Rotate section, click Custom.
  - Red arrows appear over the gel image.
- 2. Rotate the red arrows to any degree between 0 and 360 by dragging them.
- 3. Right-click the gel image and click Rotate to set your gel in the new position. Click Cancel if you do not want to set the rotation.

# Cropping a Gel Image

You can save crop settings and use them to crop other images. This feature is useful when you want to crop the same area in several images.

#### To crop a gel image

- Click Crop. A red box outlines the image area.
- 2. Drag the red box to surround the image area you want to keep.
- 3. (Optional) Right-click the image to open the Crop menu and click Save Crop Settings.
  - The Save Crop Settings dialog box appears.
- 4. (Optional) Type a name for the crop settings and click OK.
- Right-click and select Crop or Cancel. Selecting Crop crops the image to the area inside the red box.

#### To crop an image using saved crop settings

- Click Crop. A red box outlines the image area.
- 2. Right-click the image to open the Crop menu and click Load Crop Settings.
- 3. Select the saved crop settings that you want to use and click Load. The red box resizes and the crop specifications appear on the image.
- 4. Right-click and select Crop. The image is cropped to the area specified in the crop settings you selected.

#### To delete crop settings

- With an image open, click Crop.
- 2. Right-click the image inside the red box to open the Crop menu and click Delete Crop Settings.
- Select the crop settings in the dialog box that appears and click Delete.

# **Inverting Data**

Use this button to change the image data from positive to negative. Invert Data is used for negative stains and zymograms. Intensity values of bands must be greater than background to perform analysis on the gel. View the gel as a 3-D projection to determine if the data must be inverted.

# **Merging Images**

Use this button to merge a chemiluminescent blot image with a colorimetric image of the same blot. If you have used colorimetric prestained standards for a chemiluminescent blot, you can acquire an epi-white light image of the blot to show the standards and a chemiluminescent image to show immunodetection. These two images can then be merged into a combined image with both signals.

Note: Merging images can have an adverse effect on quantification. If accurate quantification is required, perform analysis on the original, separate images. Only images of the same size can be merged.

# Lane and Bands Tools

The Lane and Bands tools enable you to identify the lanes and bands in your images.

#### To open the Lane and Bands tool

In the Analysis Tool Box, click Lane and Bands, then select the tab for the Lanes or the Bands tool.



# **Detecting Lanes**

#### To detect the lanes in the image, do one of the following:

- Click Automatic if the gel image is fairly typical.
- Click Manual to detect a specific number of lanes or if automatic lane detection did not find all the lanes.

# **Using the All Lanes and Single Lane Tools**

#### To use the All Lanes and Single Lane tools

► Click the lane tool first, then click the lane to which you want to apply the change.

#### All Lanes Tools

You can modify all the lanes in your image using these tools.

- **Resize** resizes the lanes in your image. Click Resize and then drag the handles of the red squares to resize all lanes.
- **Adjust** adjusts the orientation of all lanes. Click Adjust and then drag a single corner of the lane frame. The Adjust All Lanes tool does not resize lane width.

You can add anchor points to the top or bottom borders of the rectangle by clicking the lane frame. Remove any unneeded anchor point by right-clicking it. By dragging these additional anchor points, you can adjust for "smiling" gels.

**Note:** You can move the entire lane frame when resizing or adjusting the lanes. Click Resize or Adjust and make the changes to your lanes. Then click anywhere in the frame and move it to the desired location.

**Tip:** On the PC, you can press the Shift key or the Ctrl key and use the arrow keys on your keypad to move the lane frame. On a Mac, press the Shift or the Command key and use the arrow keys.

**Delete** — deletes all lanes.

**Width** — changes the width of all lanes at the same time. Click Width and then drag one of the anchor points on any lane to change the width of all lanes.

**Tip:** You can use the plus key (+) on your keyboard to increase the lane width, and the minus key (-) to decrease the width.

You can also change the lane width by specifying the size of the lanes. Click Width, enter a number (in mm) in the box, and click Apply.

Note: You might see that the number in the box changes after you click Apply. For example, you enter 5.75 and the number changes to 5.71. The image is made up of pixels and the Image Lab™ software can only draw a boundary for the lane between pixels. If the number you specify would cause the boundary to fall anywhere on the pixels, Image Lab moves the boundary so that it falls between two pixels and updates the number accordingly.

Note: The All Lanes Width tool makes the width of all lanes uniform. Therefore, if you used the Single Lane tool to change the width of an individual lane, this change is overridden by the All Lanes Width tool.

### Single Lane Tools

Specific information about how to use each lane tool is described in the following section. You can modify an individual lane in your image using these tools.

**Add** — adds a lane to a gel image. Click Add, then click within the lane frame where you want to place the new lane. The lanes are automatically renumbered.

Note: To add a lane outside the frame, add a lane inside the frame and click Move to expand the lane outside the frame's boundaries.

**Delete** — deletes a lane. Click Delete, then click either the lane or its lane number. The lanes are automatically renumbered.

**Bend** — bends a lane to better fit the gel image. Click Bend, then drag square anchor points to fit the lane to the image.

Note: To add anchor points, left-click within the lane. Drag these anchor points to adjust the lane to fit the gel image. To remove an anchor point, right-click on it.

- **Move** moves a lane to a new position on a gel image. Click Move, then click the lane you want to move. Drag it to a new location. The lanes are renumbered according to their new positions.
- **Width** changes the width of a lane. Click Width, then click within the lane. Click on the anchor points to adjust the lane width.

You can also change the lane width by specifying the size of the lanes. Click the Width button, enter a number (in mm) in the box, and click Apply.

**Note:** You might notice that the number in the box changes after you click Apply. For example, you enter 5.75 and the number changes to 5.71. The image is composed of pixels and the software can draw a boundary only in the lane between pixels. If the number you specify would cause the boundary to fall anywhere on the pixels, Image Lab moves the boundary so that it falls between two pixels and updates the number accordingly.

# Lane Background Subtraction

#### To perform lane-based background subtraction

Select Enable Subtraction in the Background Subtraction field. Use the Lane Profile view to see the subtracted lane background.

**Disk Size** — specifies the size of a hypothetical rolling disk (between 0.5 and 99.5 mm in 0.5 mm increments) that removes background levels along the length of the lane. The size of the disk determines how closely the background level follows the intensity profile.

A large disk follows the profile trace less closely, touching fewer points along the trace and removing less background. A smaller disk more closely follows the profile trace, removing more background.

A disk radius that is too large will result in poor background removal. A disk radius that is too small might subtract actual data. For most samples, a size of ≤10 mm is usually appropriate. You can perform this task several times until you are satisfied with the amount of background removed. Use the Lane Profile tool to evaluate the appropriate disk size for background subtraction.

**Apply to selected Lane** — applies the specified level of background subtraction only to the selected lane. This option enables you to set different background subtraction levels for each lane.

# **Copying Lanes**

You can copy the lanes from one image into any other image. The entire frame and all the lanes are copied. Individual lanes cannot be copied.

#### To copy lanes between channels

- Open the Lane and Bands tool.
- 2. Select the channel that contains the lanes you want to copy.
- 3. From the Edit menu, select Copy.

The message Copying All Lanes appears.

- 4. Select the channel that you want to copy the lanes into.
- 5. From the Edit menu, select Paste.

A channel can contain only one lane frame at a time. Therefore, if you paste lanes into a channel that already contains a lane frame, you are prompted to confirm the deletion of the existing lanes. If you click Yes, the existing lanes are deleted and replaced with the copied lanes.

After the lanes are copied into the channel, you can manipulate individual lanes using the lane tools so that they are correctly positioned.

# **Detecting Bands**

Bands are detected for individual images.



#### To detect bands in the image


Click Detect Bands to open the Band Detection dialog box. Select band detection sensitivity and the lanes to which it applies.



#### **Displaying Advanced Detection Options**

You can also select Display Advanced Options to set specific parameters for band detection. When initially expanded, the values in Advanced Options are set in relation to the band detection sensitivity that you select in Detection Settings. You can set specific parameters for the sensitivity level and apply the parameters to all lanes or to a specific lane.

If later you change the band detection sensitivity in Detection Settings, the values in Advanced Options change in relation to the new sensitivity levels. You might need to reset these parameters to the original values.

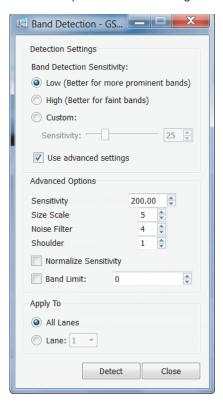

#### **Advanced Detection Options**

**Sensitivity** — determines the minimum optical density that will be defined as a band. The higher the sensitivity value, the lower the minimum signal intensity, therefore the more bands will be detected.

If the sensitivity is set too high, background staining might be detected as bands. If the setting is too low, bands of interest might not be detected.

The default sensitivity setting is 10.0. If the gel has faint bands (for example, if the optical density is less than 0.05, and counts are less than 2,000), you may want to increase this value to 20.0.

**Size Scale** — distinguishes between trends in signal intensity and random intensity fluctuations. It is the number of pixels in a vertical column that are taken together to determine whether a band is present.

The Size Scale parameter uses the size of objects in the image to determine the nature of those objects. If a gel image has high levels of background noise, a larger size scale is appropriate. At low noise levels, a smaller value is preferable. You can also increase the size scale if the gel has only a small number of thick bands scanned at high resolution.

**Noise Filter** — minimizes the number of small fluctuations (or noise) in the image that are called bands while still recognizing larger features (the real bands). This filter becomes especially important at higher sensitivity levels.

The noise filter value refers to the size of the filter in pixels (for example, a value of 2.50 equals a filter size of 2.50 x 2.50 pixels). Features smaller than the filter size will not be recognized as bands. Entering a noise filter size of zero turns it off completely.

If band detection identifies doublets as single bands, decrease the noise filter setting and/or increase the sensitivity level.

Tip: You can also decrease the Size Scale parameter instead of the noise filter to improve the detection of closely spaced bands. However, if you decrease both the noise filter and the size scale, the fuzziness around bands might be mistakenly detected as separate bands.

**Shoulder** — band detection tries to distinguish shoulders as separate bands. When looking at a lane trace, these bands appear as flat or gently sloping abutments to darker, better-defined bands (that is, there is no dip on the trace between the two bands). Increasing the shoulder sensitivity results in more shoulders being detected as bands. Changing this setting to zero results in no shoulders being recognized as separate bands.

If band detection calls a doublet a single band, check the lane trace to see if there is a dip between the peaks of the two bands. If there is no dip, increasing the shoulder sensitivity value will help resolve the two bands.

**Normalize Sensitivity** — compensates for differences in intensity between lanes.

The intensity of each lane is determined by the darkest band in that lane. For example, suppose that in all but one of the lanes the darkest band has an intensity of 50,000 counts. In the light lane, the darkest band is only 25,000 counts. With normalization, band detection will be twice as sensitive when processing the light lane, improving the detection of faint bands.

**Note:** It does not normalize for band quantitation.

**Band Limit** — enables you to limit the number of bands that will be detected in each lane, thus reducing the need for later editing.

# **Editing the Detected Bands**

You can optimize the bands in your images using these tools.

**Add** — manually add a band to a lane. Click Add, then click at the desired location in a lane. Image Lab then locates a faint band close to where you clicked.

**Tip:** You can darken your entire image to view faint bands more easily using the sliders in the Image Transform dialog box. For instructions, see Image Transform on page 92.

#### 6 | Analyzing Images

- **Delete** deletes a band from a lane. Use this feature to remove bands that are not relevant to your analysis. Click Delete, and then click on the band you want to remove.
- Adjust adjusts the height of a band. Click Adjust. Boundary lines appear on either side of each band. Move the cursor over a boundary line until you see a double-headed arrow. Move the boundary line up or down. The center is recalculated and the band appears there.


Note: You can also adjust band boundaries in the Lane Profile view.

# Molecular Weight (MW) Analysis Tools

Molecular Weight Analysis tools enable you to determine molecular weight (or base pairs, if using nucleic acid gels) by comparing a test sample with known standards.

**Note:** Before you can use the Molecular Weight Analysis tools, you must detect the lanes and bands in your image.

You can view each band's molecular weight in the molecular weight column of the Lane and Band tab in the Analysis Table.

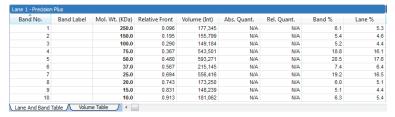

You can also display the molecular weight of the bands on the gel image by opening the Display Gel Options window and selecting Mol. Wt. from the dropdown list in the Band Attributes section. (See Displaying Gel Images on page 90 for information about displaying band attributes.)


# **Molecular Weight Standard**

You can change the standards to ones that are relevant to your samples.



#### To change standards

Click Change to access the Manage Standards dialog box. Choose another standard or add your own standards.

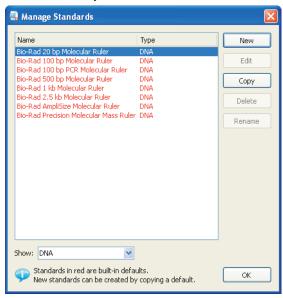


#### Standard Lanes

Standard samples are placed in the first and last lanes by default. You can specify other standard lanes by selecting the box below each lane or by entering the standard lane numbers separated by commas in the MW Analysis Tools dialog box. In the Lane and Bands view, Std appears below the lanes, identifying them as standard lanes. In the Molecular Weight Analysis view, these lanes are indicated with a check mark below the lane.

There are three standard lanes in the screen shot below (lanes 1, 16, and 17). Standard lanes are identified with a check mark in the boxes below the lane. The molecular weights of the bands in the standard lanes appear on either side of the lane frame. The red lines running from one end of the lane frame to the other identify the location of the bands in the standard lanes. You can use these lines to see where the bands in the other lanes fall relative to the bands in the standard lanes.



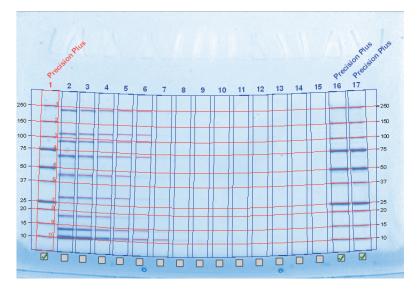

#### 6 | Analyzing Images

The values for the molecular weight in any standard lane appear in bold. In this example, the values for the first lane are in bold. (Lanes 16 and 17 are not included in the screen shot, but their values would also be in bold.) The molecular weight of the bands in the remaining lanes is calculated relative to these standards.

Lane 1 - Precision Plus		Lane 2		Lane 3		Lane 4	
Band No.	Mol. Wt. (KDa)	Band No.	Mol. Wt. (KDa)	Band No.	Mol. Wt. (KDa)	Band No.	Mol. Wt. (KDa)
1	250.0	1	203.8	1	207.9	1	210.9
2	150.0	2	118.0	2	119.7	2	121.6
3	100.0	3	99.0	3	100.4	3	102.0
4	75.0	4	70.6	4	71.6	4	72.4
5	50.0	5	45.1	5	45.8	5	46.3
6	37.0	6	26.8	6	27.2	6	27.5
7	25.0	7	17.4	7	17.6	7	17.8
8	20.0	8	12.9	8	13.0	8	13.2
9	15.0	9	10.6	9	10.7	9	10.8
10	10.0						

### **Regression Methods**

There are four regression methods.

- Linear (semilog)
- Point-to-point (semilog)
- Logistic
- Cubic spline

See Regression Calculation Methods on page 203.

# **Quantity Tools**

You can quantify bands in test samples automatically using either the Relative or Absolute tab under Quantity Tools.

#### **Relative Quantity Tab**

To compare the relative quantities of bands


1. Select the Relative tab.



2. Click Select.

#### 6 | Analyzing Images

3. Click the band you want to use as a reference. A small R appears near the band you selected.

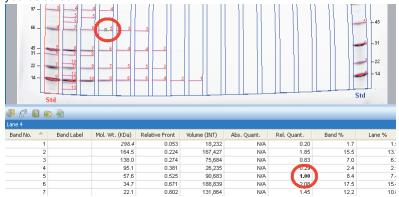

#### To review the relative quantities of bands

Go to the Rel. Quant. column of the Analysis table (Lane and Band tab). The relative quantity is the ratio of the band volume divided by the reference band volume.

All other bands now display numerical values that are relative to the reference band. If the reference band value is 1.00, values higher than 1.00 indicate that the band quantity is greater than the reference band. Values lower than 1.00 indicate the band quantity is less than that of the reference band.

# **Absolute Quantity Tab**


Absolute quantification is used to quantify bands based on known standard bands using a calibration curve.

### To calculate the absolute quantities of the bands

Select the Absolute tab.




- Click Select.
- 3. Select at least two standard (known) bands and assign quantity values. Small A's appear (shown below circled in red) near the bands you selected.

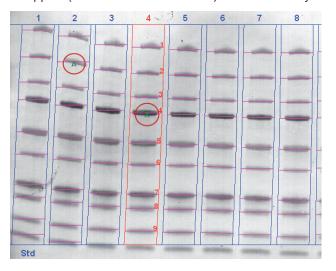

#### 6 | Analyzing Images

The values display in the Standard Bands table. The greater the number of known bands and the wider the range of their values, the more accurate the absolute quantity calculation of the unknown bands will be.

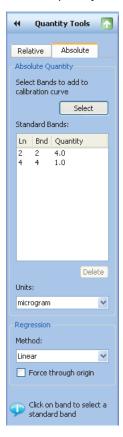

Note: Any standard band selection can be deleted. To do so, select the entry in the Standard Bands field and then click Delete.

- Select a unit of measure from the Units dropdown list.
- 5. Select a regression method from the dropdown list.

Consider the following guidelines when making your choice.

**Linear** — generates a straight line that is the best fit of the values you provided and is preferred in most cases.

**Point-to-point** — generates a curve in which each data point is connected directly to the next, regardless of the shape of the resulting curve.

**Cubic spline** — generates a smooth curve that connects each data point. At least four standard points are required to use this method of least-squares polynomial fits.

**Table 7. Regression Methods** 


Regression Method	Minimum Number of Standard Bands	Minimum Number with Force Through Origin Option
Linear	2	1
Point-to-point	2	1
Cubic spline	5	4

6. Click Standard Curve on the toolbar and select the Absolute Quantity Standard Curve tab.

#### 6 | Analyzing Images

The calibration curve displays. Standards are represented by green triangles. Unknown values are represented by red triangles.

Note: Selecting the Force Through Origin checkbox always starts the standard curve graph at 0,0, regardless of the best curve fit.

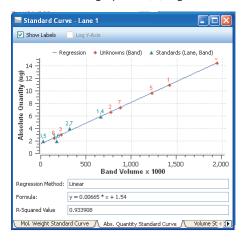

Note: Clicking the Standard Curve table generates a crosshair tool that displays the numerical values associated with the placement of the cursor in the graph.

# **Annotation Tools**

You can annotate results with text and arrows to emphasize areas of interest.



#### **Add Annotations**

**Text** — enables you to add text annotations to images to emphasize important details. Click Text, then click an area on the image where you want to insert your comment. A box appears with a dotted-line border. Type your comment in the box.

**Tip:** To add a new line to the box, place your cursor where you want the break and press Shift+Enter.

#### To move the box

- Click and drag the box to change its position.
- **Arrow** enables you to add arrows to images to emphasize important details. Click where you want the arrow to start and drag to stretch the arrow point to the location you want to emphasize. To move the arrow on the image, click the middle of the arrow and drag it to the new position.

#### To change where the arrow points

Click either end of the arrow. Square boxes appear. Drag a box to change the length or orientation of the arrow.

# **Alignment**

The alignment buttons enable you to align multiple annotations, such as lane numbers, which you have manually added.

#### To select multiple annotations

Press the Ctrl key (Command key on a Mac) and click each item or drag a selection box around them.

# **Text Properties**

You can change the type size and font of your text annotations.

- **Font** click the box you want to change. Open the dropdown Font menu to show all fonts installed on your system. Select a new font for the text annotation.
- **Size** click the box whose font you want to resize. Open the dropdown Size list to increase or decrease the size of the text. You can set the font size from 6 to 72 points using the dropdown list.

# Color

You can change the color of text annotations to make them visible with any color scheme and emphasize them further by adding a color to the annotation's background, which is invisible by default.

#### To change the color of multiple items

- Press the Ctrl key and click each item.
- **Foreground** click a text annotation or arrow. This activates the Foreground field, so you can select a foreground color from the dropdown list.
- **Background** click a text annotation. This also activates the Background field so you can select a background color from the dropdown list.

### Rotate

You can rotate text annotations 90° to the left or right by clicking the Rotate buttons.

# **Volume Tools**



Volume tools enable you to manually quantify features on a sample image when automated lane and band analysis is not appropriate or possible, such as in dot blots.



You can use Volume tools to quantify the signal intensity of bands, spots, arrays, and other image data. Define an area of interest by surrounding it with a shape. You can choose a rectangle, circle, freehand, or lane shape by clicking the appropriate button under the Volumes field.

A default label appears within the shape drawn. The volume label is assigned a sequential number and can be one of three types:

- U unknown
- Std standard
- B background



Each new volume you create initially has a red border, which indicates that the volume is selected. When you click elsewhere on the image, the border changes to blue, indicating that the volume is no longer selected.

**Note:** Double-click a volume area to change its properties.

#### To review data for the volumes

Open the analysis table and select the Volume tab. Volumes are listed based on their number and/or the associated information per volume. See Volume Measurement Definitions on page 100.

#### **Volume Types**

You can define the volume type (unknown, standard, or background), the quantity of standard volumes, or enter a custom name to replace the default label.




Unknown volumes are volumes you want to quantify.

Standard volumes are used for absolute quantities. See Absolute Volume Quantity on page 140.

Background volumes are used to remove the background from the calculation. The result of volume background subtraction appears in the Adjusted Volume column of the analysis table (Volume Table tab).

Note: This volume type needs to be assigned only when using Global Background subtraction.

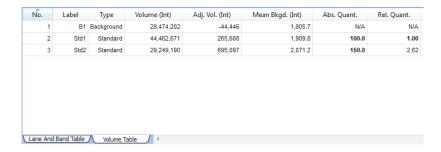

#### **Volume Background Subtraction**

When you draw a volume, some nondata background pixels might be included inside the volume. These background pixels usually have an intensity value that you do not want to include in your volume quantification. There are two methods of calculating this background intensity for your volumes: local and global.

- **Local** local background subtraction calculates a separate background intensity for each unknown and standard volume you create. For each volume, the intensities of the pixels in a 1-pixel border around the volume are added together and divided by the total number of border pixels. This gives an average intensity for the background around each, which is then subtracted from the intensity of each pixel inside the volume. If the background value is greater than the pixel value inside the volume, the background-adjusted quantity of the volume could be <0. In this case, redraw the border for this volume.
- **Global** global background subtraction calculates a single background intensity for the entire gel. This average background intensity is then subtracted from all the volumes in the gel. The average intensity of the pixels in the background volume is calculated and subtracted from each pixel in all standard and unknown volumes. Therefore, it is not necessary for the background volume area to be the same size as your unknown.

#### To calculate global background subtraction

- Use one of the Volume Tools to create a volume in a representative background region of your image (that is, a nondata region similar to the background surrounding your data).
- 2. Double-click the volume. This opens the Volume Properties dialog box.
- 3. Select the Background option button.

#### Note:

If you select Global in the Volume Tools toolbox but do not define a background volume as described, no background subtraction is performed.

- If you create more than one background volume, all the pixels in those background volumes are used to calculate the average background. Your background volume(s) will have default names B1, B2, and so on based on the sequence in which they were created.
- If the region you defined as background has a higher average intensity value than your data object, you obtain a negative value for your adjusted volume in the analysis table. If this happens, select a new background region with less intensity than your data object.

#### **Relative Volume Quantity**

You can choose any one volume as a reference volume by selecting the Reference Volume checkbox in the Volume Properties dialog box. The reference volume is indicated by an asterisk on the volume label, for example, U1\*.

Relative quantities are displayed in the Relative Quantity column in the analysis table (Volume Table tab). The relative quantity is the ratio of the backgroundadjusted volume divided by the background-adjusted reference volume.

All other volumes now display numerical values relative to your reference volume. Values higher than 1.00 indicate that the volume is greater than the reference volume. Values lower than 1.00 indicate the volume is less than the reference volume.

# **Absolute Volume Quantity**

If you have drawn your volume around an object of known quantity, you can use it to calculate the quantity of your unknown volumes. The quantities of your unknown volumes are calculated based on the standard volumes and the selected regression method.

#### To classify a particular volume as a standard

- 1. Double-click the volume. This opens the Volume Properties dialog box.
- 2. Select the Standard option button and enter the quantity in the Quantity box.
- 3. Click OK to close the dialog box.

Standard volumes will have the default names S1, S2, and so on, based on the sequence of their creation.

#### To review the regression curve

Open the Standard Curve window and select the Volume Standard Curve tab.

#### **Regression Methods**

Three regression methods are available to generate the volume quantification curve used for absolute quantity: linear, point-to-point, and cubic spline. To display the standard curve, click the Standard Curve button on the toolbar and select the Volume Standard Curve tab in the Standard Curve dialog box. See Regression Calculation Methods on page 203 to learn how each of these methods is calculated.

The data for volume standards are found in the Absolute Quantity column of the Volume Table.

Note: Selecting the Force Through Origin checkbox always starts the standard curve graph at 0,0, regardless of the best curve fit.

#### **Alignment**

Align volumes by using the appropriate alignment button. To select several volumes, Ctrl-click each one, then select one of the alignment buttons. Hover over any of the six alignment buttons to display its function (Align Left, Align Right, etc.)

Copy and paste selected volumes by pressing Ctrl+C to copy. Press Ctrl+V to paste.

When you click the Standard Curve button on the toolbar, a chart displays all unknown and standard quantities.

# **Generating Reports**


After viewing results, you can generate a report that displays the analyzed gel images, all of the settings used in the protocol, and as much information about the data as you want to include.

You can choose print settings within the Report Settings dialog box in the Edit menu or by clicking Report in the main toolbar.



#### To produce a preview of your report

Click Report on the toolbar.

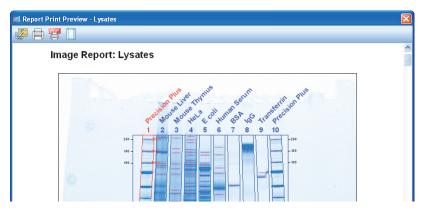


Use the following dialog boxes to customize the content in your reports. Doing so does not modify the data from the analysis.

# Report Settings

Use this dialog box to customize the content of your report.

#### **General Tab**

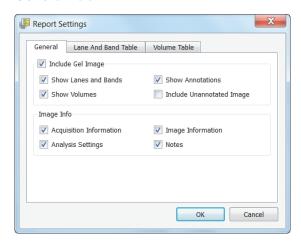

The General tab includes the following settings:

**Include Gel Image** — specify whether the image is included in the report.

If the image is included, the following options determine which overlays are displayed on the gel image:

- Show Lanes and Bands П
- **Show Volumes**
- **Show Annotations**
- Include Unannotated Image

When you select this checkbox, both the annotated image (if available) and unannotated images are printed in the report. The unannotated image precedes the annotated image in the report.


Note: If you disable the other checkboxes in the Include Gel Image group, the Include Unannotated Image checkbox is automatically disabled.

	Image Info —	specify what	information i	is included	in the report.
--	--------------	--------------	---------------	-------------	----------------

- Acquisition Information
- Analysis Settings
- Image Information
- Notes
- **Signature History** details on when and why a secure document was signed

If a secure document has been signed, the user name, the date and time of the signature, and the reason for signing are all included in the Signature History section. If a document has not been signed, this section is omitted from the image report.


### Lane and Band Table Tab

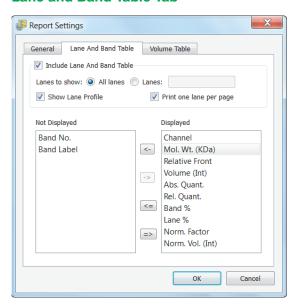

The Lane and Band Table tab includes the following settings:

- **Include Lane and Band Table** specify whether to include the Lane and Band table in the report.
- **Lanes to show** specify which lanes to display in the report.
- **Show Lane Profile** specify whether to display the Lane Profile view.
- **Print one lane per page** specify whether all lanes are printed on one page or whether each lane is printed on a separate page (adds a page break after each lane).
- **Show Lane Profile** include the lane profile for each lane.
- Not Displayed/Displayed remove columns that you do not want to display in the report. By default, the report displays all columns from the Lane and Bands table.

### **Volume Table Tab**

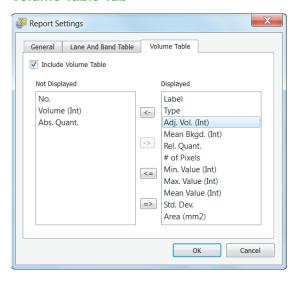

The Volume Table tab includes the following settings:

- **Include Volume Table** clear to exclude this information from your report.
- Not Displayed/Displayed remove columns that you do not want to display in the report. By default, the report displays all columns from the Volume table.

# Print Report

Click the Print Report button to print your report.

# Print Report to a PDF File

The Print Report to .pdf File button opens a Save dialog box so the PDF file can be saved on your system.

# Adjust the Printer Settings

The Printer Settings button accesses options for paper size, orientation, and page margins.

### **Exporting Results** 8

The most convenient way to archive complete information about experiments is to produce reports. However, you might want to export only gel images or analysis table data for analysis in different programs, such as Quantity One®, FPQuest™, or InfoQuest™FP software. Or you might need exported files for presentation or publication.


# **Exporting Gel Images**

Image Lab™ software includes features for exporting gel images several ways.

- You can export displayed image data to a publication (choose Export for Publication).
- You can export raw image data as a 16-bit .tif file (choose Export for Analysis).
- You can export image data to PulseNet. Doing so reduces the image to an 8-bit .tif file, limits its resolution, and restricts its file size to 300 Kb.
- You can export lane and band tables as well as volume tables to a spreadsheet program or to a file.

#### 8 | Exporting Results

The options to export gel images are available using the Export option in the File Menu.

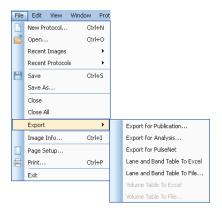

# **Exporting Gel Images for Publication**

Use this format only to export visual information to presentation or word processing software, such as PowerPoint or Word.

### To export a displayed image to a file

Select File > Export > Export for Publication.


Note: You can select from .bmp, .png, .jpg, and .tif formats. The gel displays with any lanes, bands, and annotations that appear on the screen.



### In this dialog box you can:

- Select the entire image or the current view
- Select the resolution or specify a custom resolution
- Specify the publishing dimensions
- View the resulting published image size and dimensions

You can zoom in on an area in a current view and export only that area, or you can export the entire image. You can exclude annotations or overlays by clicking Display Gel Options on the toolbar to access the appropriate settings.

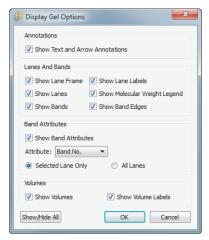

# **Exporting Gel Images for Analysis**

### To export an image for analysis

Select File > Export > Export for Analysis.

This exports the raw data only as a 16-bit .tif file.

**Note:** 16-bit .tif images are not compatible with all image viewers.

The image might require contrast adjustment when it is imported into analysis software. This option creates a file that can be analyzed in other programs such as Quantity One, FPQuest, or InfoQuestFP software.

# **Exporting Gel Images to PulseNet International**

### To export an image to PulseNet International

Select File > Export > Export for PulseNet.

Image Lab reduces the image to an 8-bit .tif image file. Resolution is limited and file size is restricted to 300 Kb.

# **Exporting Lane and Band Tables to Excel**

If you have Excel (or Numbers on a Mac) installed on your computer, you can export the data to the spreadsheet application.

### To export the data to Excel (or Numbers)

Select File > Export > Lane and Band Table to Excel.

This opens a table directly in the spreadsheet program. You can then use the Save As option to produce other formats.

# **Exporting Volume Tables to File**

### To export an image as a CSV file

Select Export > Volume Table to File.

Image Lab exports the image as a comma-separated values (CSV) file so the data file can be opened in a database application.

## Screenshot Tool Export

Use the Screenshot tool, available on the toolbar, to capture a displayed image to the clipboard or to save it to a file (.bmp, .gif, .jpg, or .png).

## Analysis Table Export

You can export table analysis data from the File menu or by using the export buttons at the top of the Analysis Table window.

The Analysis Table window has several buttons to export data to different formats, depending on how the data are to be presented.

# Copy Analysis Table to the Clipboard

Copies the analysis table to the clipboard and then pastes the analysis table into word processing or presentation applications. It is best to use the vertical table orientation when copying to an 8.5 x 11-inch page to accommodate the columns.

# **Export Analysis Table to a File**

Exports an analysis table as a CSV file, so your data file can be opened in a database application.

# **Export Analysis Table to a Spreadsheet**

Enables you to use sort and formula functions of a spreadsheet program with your data. If you have Excel (or Numbers on a Mac) installed on your computer, the data open in the spreadsheet program.

# **Image Lab Logs**

# **Image Lab Logs**

Image Lab™ software provides three types of logs.

- **Instrument log** records events related to the instrument, including calibrating the instrument and the success or failure of the calibration. This log file is visible only if the computer running Image Lab is connected to an instrument.
- **System log** records events related to running Image Lab, including enabling or disabling of secure mode and the users who log on to or log off of Image Lab.
- **Document log** (Security Edition only) records events related to the creating and modifying of secure protocol and image files.

# **Viewing the Instrument Log**

Instrument operations that you perform with Image Lab are recorded in the instrument log.

### To open the Instrument Log Viewer

From the View menu, click the log for your instrument.



The menu displays the specific instrument to which you are connected. The GS-900™ densitometer is shown as an example, above.

Note: You can view an instrument log only if Image Lab is connected to the specified instrument.

The instrument log includes the following information.

- date and time of the action
- the user name of the user logged into the computer at the time of the action
- the user's security level
- the domain where the current user is logged in
- the type of action performed
- the result of the action taken, including complete records of the calibration
- the user's reason for the action, if provided




# Viewing the System Log

Events related to running Image Lab, including enabling or disabling of secure mode and users logging on to or logging off from Image Lab, are recorded in the system log.

### To open the System Log Viewer

From the View menu, click View System Log.

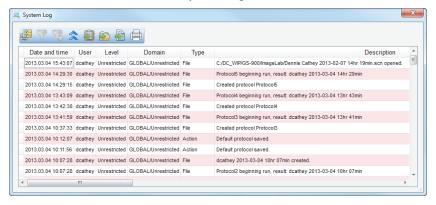

# **Viewing the Document Log**

The View menu displays a list of logs for each open document. The document log captures information about creating and editing the Image Lab protocol and image files.


Note: Document logs are viewable only in Image Lab Security Edition, and only on Windows-based computers. Document logs are not viewable on the Mac.

The document log captures changes from:

- Image tools
- Lane and bands tools
- Normalization tools
- Molecular weight tools
- Quantity tools
- Volume tools


You can view the document log for any open file. This file can be a previously saved file, or it can be a newly created protocol or image file that is open on your desktop but which has not yet been saved.

If the document was previously saved, the log is identified with this name. If the document has not been saved, the log is identified with the time stamp, showing when the document was created.

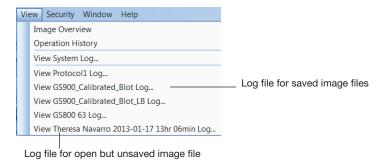

### To open the Document Log Viewer

From the View menu, click the document log that you want to view. A typical document log is shown below.

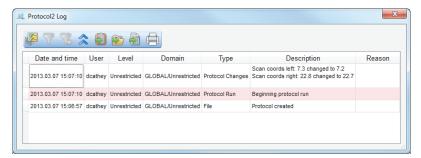

# **Displaying Log Data**



The display toolbar is located above the log.

# Displaying Data Columns in Logs


By default, the logs display the following columns:

- Date and time
- User
- Level the security role of the user
- Domain domain where the current user is logged in
- Type the type of event
- Description the actual event captured
- Reason (For Secure Mode only) the reason for signing a document

You can change your view of any log by displaying or hiding data columns.

### To display or hide columns

- From the View menu, open the log file.
- 2. Click the Display log viewer options icon.
- 3. In the Display Column Options dialog box, use the arrow keys to move columns between the Not Displayed and Displayed lists.

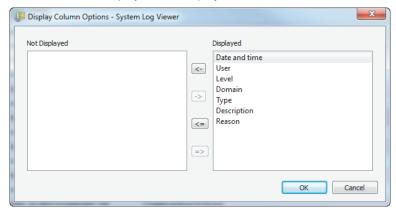

Click OK.

# Filtering Data in Logs 🔀 🍞



For all Image Lab logs, you can filter the entries in the following columns:

- Date and time
- User
- Type

For example, you can set the filter in the Type column to File and Image Lab will display only the rows where Type is equal to File.

### Setting filters in logs


### To filter the entries of a column

In the log, click the heading of the column that you want to filter. The Display Filter options icon in the Actions tool bar is enabled. (By default, the log opens with all event types displayed.)

Note: The Display Filter options icon is enabled only when you click on a column that can be filtered. The Display Filter options icon remains disabled if you click a column that cannot be filtered.

2. Click the Display Filter options icon to display the Select filter values dialog box.

**Tip:** Alternatively, right-click the heading of a filterable column to display the Select filter values dialog box.

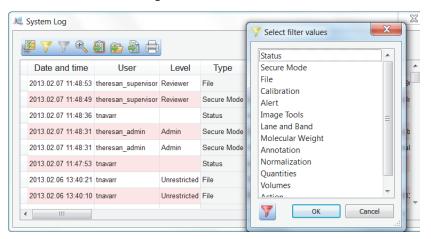


Select a value and click OK.

**Tip:** You can select multiple values to filter.

### Removing filters in logs

The Remove all filters icon is enabled only after you set a filter on a column. This icon clears all filters on all columns.

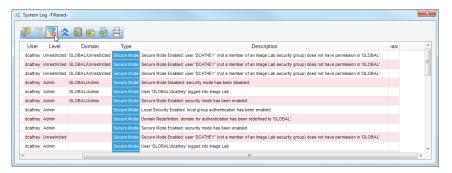

Each filterable column has a Remove filters icon as well. This icon removes the filter for that column only.



# Collapsing or Expanding Data Rows 🚨 🎽





You can expand the size of rows in any log to display the full content of the row, or you can collapse the row size to display more rows in the table. The icon toggles between collapse and expand.

### To display the full content of a row

► Click the Collapse or expand rows icon to expand the row height. The row adjusts to fit the text of the longest entry.

This action expands all rows in the log where an entry is longer than the width of the column. Rows that do not need to wrap are not affected.

### To display more rows in the table

Click the Collapse or expand rows icon to collapse the row height to the default height setting.

This action collapses all rows in the log to display data on a single line.

# **Exporting Logs** 🚨 🔁 🖺



From the log viewer, you can:


Copy log entries to the clipboard — copies the log entries to the clipboard, enabling you to paste them into a word processing or presentation application.

**Tip:** You can also open a log in the log viewer and press Ctrl+C to copy the log entries to the clipboard.

- **Export log entries to a file** exports the log entries as a CSV file that can be opened in a database application.
- **Export log entries to Excel** exports the log entries to an Excel file where you can use Excel's sorting and formula functions to manipulate your data. If Excel is not installed on your computer, this feature is disabled.

### To export a log file

- From the View menu, click the log you want to view and export.
- 2. Click one of the icons to export the data in one of the supported formats.

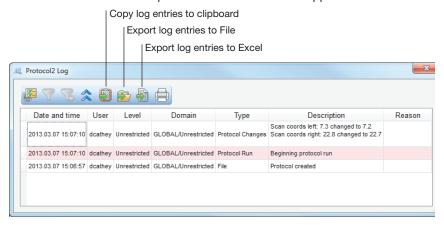


# Printing Logs

Log data can be sent to a printer or saved to a PDF file.

### To print a log

- From the View menu, click the log you want to view and print.
- 2. In the log viewer, click Print log.

The Log Print Preview window displays the contents of the log file.

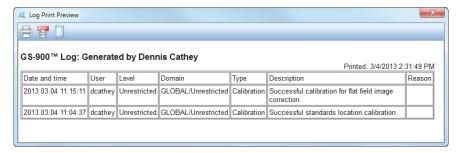

From the Log Print Preview window, you can:

- Click Print log to print the log to a printer.
- Click Print log to PDF to save the log to a PDF file.
- Click Adjust printer settings to prepare the file for printing.

# 10 Using the Security Edition

### 21 CFR Part 11

Image Lab™ Security Edition is a module within the Bio-Rad Image Lab software that assists users in meeting the Food and Drug Administration's regulations on good lab practices in the pharmaceutical and biotechnology industries. The Security Edition enables system administrators to ensure that Image Lab operates in compliance with Title 21 of the Code of Federal Regulations (CFR) Part 11 within a closed system. A closed system is defined as "an environment in which system access is controlled by the persons who are responsible for the content of electronic records that are on the system" (Section 11.3 (b) (4).

### Note:


- Image Lab Security Edition is not supported by the Mac.
- The security controls built into Image Lab Security Edition must be properly configured and administered by the system administrator(s) in your organization in order to be secure and in compliance with 21 CFR Part 11.
- Bio-Rad makes no claim that Image Lab Security Edition software is CFR-compliant in and of itself, nor does the company guarantee compliance for the user. Your organization must establish policies and standard operating procedures that work in conjunction with the tools provided by Bio-Rad to ensure compliance with 21 CFR Part 11.

## Standard Mode versus Secure Mode

Image Lab Security Edition can be run in either of the following modes:

- **Standard Mode** in this mode, there are no restrictions on controlling the instrument, operating the software, or changing the documents.
- **Secure Mode** in this mode, in addition to the features available in standard mode, extensive security features are enabled, including document signing and log creation.

When Image Lab is in secure mode, a padlock symbol appears in the left corner of the status bar. If no padlock is present, the software is running in standard mode.

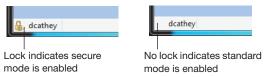

When Image Lab is installed, by default it is set to run in standard mode. It continues to run in this mode until a user with Image Lab Administrator privileges enables secure mode.

This chapter assumes you are running the application in secure mode unless otherwise noted.

# **User Names, Groups, and Roles**

When Image Lab Security Edition is run in secure mode, you must log in with a user name and password in order to run the application. The Microsoft Windows system administrator generally creates the user names and passwords. The Image Lab administrator defines the groups to which the Security Edition roles will be associated.

Note: See Setting Up Users and Groups on page 183 for additional information about setting up groups, user names, and passwords.

There are four default roles in Image Lab Security Edition. Each role is associated with one of the four default Image Lab user groups, and each user is assigned a role that provides the user access to specific features in the software. Table 8 lists the default Image Lab Security Edition user groups and the corresponding roles. It also provides a brief description of the permissions for each role.

**Note:** System administrators can change the user group names, if necessary, to meet their local company standards.

Table 8. Image Lab Security Edition groups and roles

User Group	Image Lab Role	Description
TDS_Administrator	Administrator	Users with this role can enable or disable secure mode. Administrator users can also view log files, but they do not have access to any other features.
TDS_User	Supervisor	Users with this role have full access to all features and functions of the application. They can also sign files. Supervisors cannot enable or disable secure mode.
TDS_Tech	Clinician	Users with this role can perform instrument operations, run existing protocols, and view protocol files, results files, and log files. They can also sign files. All other access is restricted. For example, they cannot create or edit protocols.
TDS_Guest	Reviewer	Users with this role can open and view protocol files, results files, and log files. They can also sign files. Users with this role cannot create new protocols and cannot change protocols. All other access is restricted.

### **Role Restrictions**

Your role determines which features of the security edition you have permission to use. If you attempt to perform an action that is not permitted for a user in your role, you will see an error message. In some instances the user's role determines which Security Edition features are visible and/or enabled. Therefore, you might not see all of the features described in this chapter.

Table 9 lists the Image Lab Security Edition functions that each role has permission to perform.

Table 9. User access to features by role

Function	Administrator	Supervisor	Clinician	Reviewer
Enable/disable secure mode	Х			
View log files	Х	Χ	Χ	Χ
Set up and recalibrate the instrument		Х		
Create new protocols		Χ		
Open existing protocols		Χ	Х	Χ
Run protocols		Х	Х	
Edit protocols		Χ		
Create new images		Χ	Х	
Open existing images		Χ	Х	Χ
Edit images		Χ	Χ	
Analyze images		Χ	Χ	
Sign documents		Χ	Χ	Χ

# **Starting Image Lab Security Edition**

When secure mode is enabled, you are prompted to log on when you start the Image Lab application.

### To start Image Lab in secure mode

- 1. Click the Image Lab icon to start the application.
- In the Log on to Image Lab dialog box, enter your user name and password.



Click OK. 3.

Note: If you have any questions or problems logging on, see your system administrator.

# **Electronic Records**

Image Lab Security Edition enables you to create secure electronic records as defined by 21 CFR Part 11. In Image Lab, the following are electronic records:

- Protocol files
- Image files
- Document log files
- Instrument log files
- System log files

### **Unsecured Documents**

Protocol and image files created in standard mode or created with previous versions of Image Lab are unsigned, unsecure documents. Unsecure documents remain unsecure. When you open these files in secure mode, you can make changes to these files and save them without restrictions. A document log is generated for changes to these files, but the log is not viewable in standard mode. You must log in to secure mode to see the log.

You can also open unsecure documents in secure mode and sign them. You can make changes to these files and save them as secure documents. The original unsecure document remains unsigned and unsecure. The new document is saved as read-only with an incremental revision number and Image Lab generates an audit log. Image Lab will not be able to overwrite the secure file, but it can overwrite the original unsecure file.

### Secure Documents

Protocol and image files created in secure mode are known as secure documents and generate a log. However, until they are signed, such documents save as standard files. Only after they are signed are they saved with the secure extension(.sptl or .sscn). Image files generated from signed protocols automatically open the signing dialog box when the scan completes.

If you create a default protocol based on an existing signed protocol, the resulting new default protocol cannot be edited.

Note: A secure document can be saved unsigned. Likewise, a signed document can be saved as a new, unsigned file. In this case, the saved file is still a secure document. All changes to the document are captured in the document log.

Documents created in secure mode can be signed at any time. When a secure document is signed, it is saved as read-only. Image Lab cannot overwrite the document. You can open signed documents and sign them again. The newly signed file is saved as a second revision, and the new signatures are captured in the document log.

**Note:** Image Lab can never overwrite any signed file.

You can open secure documents in standard mode of Image Lab. Selecting Save As from the File menu creates an unsecure file. The original file is preserved. No log is generated for the new, unsecure document. However, an entry is added to the original log that the file has been reverted to unsecure.

#### Secure Documents in SAM mode

Images saved in signal accumulation mode (SAM) follow the same rules as the main image. The rules apply when the image is saved from within a protocol. If you create an image from a signed protocol, all resultant images must also be signed. If you try to save a file more than once, a warning message appears asking you to confirm the re-save. And if you confirm this re-save and click OK, Image Lab creates a new file with an incremental revision number.

# **Modifying Secure Documents**

You can open and change a signed document in Image Lab. The original document is read-only and cannot be overwritten. Saving the revisions opens the Save As dialog box. The changes are saved into a new revision with an incremental revision number.

Each time a secure document is modified, you must provide a reason for each change before you can sign the document. The modifications are logged in the document log. The new signed document takes with it the entire history of the original document in its log.

# **Signing Documents**

### To sign a document

- Select the protocol or results file.
- 2. From the Security menu, click Sign Document. The Signing Document dialog box appears.



Enter the user name and password of a user authorized to sign documents.

**Note:** The user name and password can be for a user other than the current user.

4. Enter a reason for signing the document. Typical reasons include review, approval, responsibility, or authorship.

Note: You must provide a reason in order to sign the document.

The user name, date and time of the signature, and reason for signing are always included in the Signature History section of the image report (see page 145).

Click OK.

A Save File dialog box appears.

Enter a new name for the file and click Save.

Note: Signed protocols are saved with an .sptl extension, signed images are saved with an .sscn extension.

# **Document Logs**

Any changes made or actions performed on a protocol or image file generate a document log documenting each change or action. This log is created as soon as the protocol or scan file is created. Protocols are updated when the protocol is saved or run. If you make changes to a file and save it as a new file, regardless of whether you are in secure mode, the document log is preserved in the new file. If you are in secure mode, the signature of the previous file is noted as part of the document log.

Note: When a protocol is run, its log entries are copied to the resulting image file.

All major actions and changes are audited (generate a document log). Examples of auditable actions include:

- Signing a file
- Changing protocol settings
- Changing the image, for example, by cropping or rotating it
- Modifying the analysis, for example, changing the lanes and bands, or adding or editing annotation to the images

Minor changes that affect only the display are not audited, such as:


- Selecting different columns in the Analysis Table
- Changing the display options in Lane Profile
- Annotating or labeling the image using Annotation Tools

Each change you make to a signed protocol or image file must be documented in the Reason for Change dialog box.

# **Viewing the Document Log**

The View menu displays a list of logs for each open document. The document log captures information about creating and editing Image Lab protocol and image files. You can view the document log for any open file. This file can be a previously saved file, or it can be a newly created protocol or image file that is open on your desktop but which has not yet been saved.

If the document was previously saved, the log is identified with this name. If the document has not been saved, the log is identified with the time stamp of when the document was created.

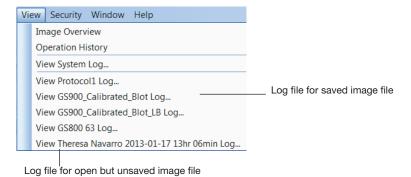

The document log includes:

- Date and time
- User
- Level the security role of the user
- Type the type of event
- Description the actual event captured
- Reason the reason for signing a document

For more information about logs, see Chapter 9, Image Lab Logs on page 155.

# A Maintenance and **Specifications**

# Cleaning the Sample Trays

Clean the sample trays with standard laboratory detergents or mild solvents such as EtOH or MeOH. Use lint-free tissue to wipe them dry. Dust particles or lint on a sample tray can glow under UV illumination.

# **UV-B Fluorescent Lamp Replacement**

The UV-B lamps provide service under normal usage for 4-5 years. In case of a lamp failure, one of the following error messages appears:

- Fault in UV bulbs Bank 1&2
- Fault in UV bulbs Bank 1
- Fault in UV bulbs Bank 2

#### To replace the lamps

- Turn off the instrument and unplug the AC cord.
- 2. Remove and save the seven screws around the outside edge of the imager back. Do not remove the screw at the bottom left outside corner of the imager back.
- 3. Carefully slide the top cover toward the back of the instrument until it is all the way off.
- 4. On the left side of the instrument a single screw holds the lamp assembly in place. Remove and save this screw.

### A | Maintenance and Specifications

Carefully slide out the lamp assembly. Although only one lamp bulb might be out, it is recommended that all lamps be replaced to ensure even illumination and to reduce the need to access the lamps again.



- Replace all lamps with UV-B fluorescent lamps (Bio-Rad part catalog #900-0217). Make sure each lamp is properly seated in its holder.
- 7. After replacing all the lamps, carefully slide the lamp assembly back into the instrument. Press on the side of the lamp assembly to the right of the screw hold-down tab, to make sure it is fully seated into the instrument. If the lamp assembly is not fully seated, the following error message appears the next time it communicates with Image Lab software: Light tray not detected.

# **Error Messages**

### Sample tray not detected

The sample tray has a magnet that senses when the sample tray is inserted. If you attempt to image and this message appears, the sample tray might not be pushed in all the way. Press the tray in until the magnet grabs the tray.

### Light tray not detected

This error message appears if the lamp assembly is not fully seated into the imager. Reseat the lamp assembly as described on the previous page.

### Door was opened during imaging

To prevent UV radiation exposure, the UV-B lamps do not turn on unless the front sample tray door is fully closed. This message also appears when the instrument cover has been removed.

A | Maintenance and Specifications

# **B** Troubleshooting

Problem	Possible Cause	Solution
Green LED on front panel is off	The AC power cord is not connected.	Connect an AC power cord to the imager and an appropriate wall socket.
	The power is off.	Turn the power switch on.
Front panel LEDs remain flashing	The firmware is not loading.	Call Bio-Rad Technical Support for help
Image is not visible on the monitor	The sample tray does not contain a sample.	Place a sample on the tray and run the protocol again.
Image is not bright enough	The gel is underdeveloped.	If initially stopped after 2 min, allow the gel to develop for 2–3 more min.
Red LED flashing on front panel	The sample tray not inserted all the way.	Push the sample tray into the imager until the magnet pulls the sample tray.
	The door interlock is broken	Close the door.
	A bulb has failed.	Replace all the lamps according to the instructions in UV-B Fluorescent Lamp Replacement on page 177.
	The communication has been interrupted.	Make sure imager power is on and the USB cable is connected to the PC. Restart the imager.
	The imaging button on the front panel was pressed when the software was not running.	The red light stops flashing when the software is opened. Press the imaging button again to initiate imaging.

## C Setting Up Users and **Groups**

## **Setting Up Image Lab Users and Groups**

This appendix explains how to set up users and groups to run Image Lab™ Security Edition software in secure mode.

Note: This task requires System Administrator privileges on the client computers (and possibly their domain) on which Image Lab is installed.

### **User Accounts**

To give users access to Image Lab Security Edition, you can create new Windows user accounts, add existing user accounts to the four default user groups specified in Table 8 on page 169, or rename any of the four default user groups.

- A user account can have any name, but must have a password defined for the account. See the section on Password Security, on page 194 for information about setting passwords for maximum security.
- Each user can belong to the Image Lab Administrator group and one other Image Lab user group.


For example, a user can belong to the Administrator group and the Supervisor group, but a user cannot belong to both the Clinician group and the Supervisor group.

## **User Authentication and Group Membership**

Authentication in Image Lab is made up of two processes: user authentication and group membership evaluation.

#### **User Authentication**

In the user authentication process, Image Lab matches (authenticates) a user name with permissions assigned to that user name on the authentication domain. That domain can reside on your local computer (a local domain) or on a network server (a network domain). You set the location of this domain in the Domain to be used for authentication field of the Security Preferences dialog box. This name must be the exact name of your local computer or network server. For details on how to find this name see Finding the name of your authentication domain, on page 185.

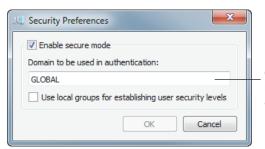

Enter the name of your local computer, or the name of the network server used for your authentication domain.

In the Security Preferences dialog box, if you (or your network administrator) choose a local domain to be used for authentication, you are considered a local user. If you or your network administrator choose a network domain, you are considered a domain user.

## **Group Membership Evaluation**

In the group membership evaluation process, Image Lab verifies that a user is a member of one or more of the four default Image Lab user groups (TDS\_Administrator, TDS\_User, TDS\_Tech, or TDS\_Guest). The valid members for each of these groups can be specified in one of two places, as defined by the Use local groups for establishing user security levels checkbox on the Security Preferences dialog box.



This checkbox specifies whether valid users and groups are defined on (1) the local Image Lab computer, or (2) on the network domain.

If this checkbox is selected, only users and groups that are defined on the local computer (on which Image Lab is installed) are recognized. If this checkbox is not selected, only users and groups that are defined on the network domain are recognized.

## Finding the name of your authentication domain

Your authentication domain can be hosted on your local computer (a local domain) or on a network server (a network domain).

#### To find the name of your local domain

- From your Start menu, open the Control Panel.
- 2. Click System.

The System window appears. In the Computer name, domain, and workgroup settings section, the name of your local computer is shown as Computer name.



### To find the name of your network domain

- From your Start menu, open the Control Panel.
- 2. Click System.


The System window appears. In the Computer name, domain, and workgroup settings section, the name of your network domain is shown as Domain.



## **Configuring Users and Groups on a Local Computer**

#### To set up users and groups on a local computer

- In the Windows Control Panel, select Administrative Tools.
- 2. Select Computer Management.
- In the Computer Management window, expand the System Tools folder, and then expand the Local Users and Groups folder.

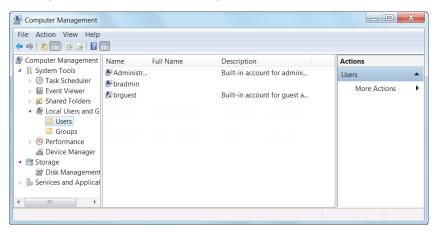

#### To create a new user on a local computer

Click on the Users folder to open it and select Action > New User. Alternatively, use the right-click context menu.

The New User dialog box appears.



- Fill out all the fields:
  - **User name** the user name must be unique.
  - **Full name** the full name must be unique.

Bio-Rad recommends using the user's actual full name, as this name will be shown in the document log and all the log reports. This is a requirement of 21 CFR 11.50a.

**Description** — this field must also be filled out.

Bio-Rad recommends entering the user's title as the description.

**Password** — enter and confirm a password for the user.


**Tip:** Select the User must change password at next logon checkbox. This prevents the Windows system administrator from knowing the passwords of the users.

Note: If you select the User must change password at next logon checkbox, the user must actually log on to Windows and change the password before using Image Lab Security Edition. Otherwise, Security Edition will not recognize the user.

#### To create a new group on a local computer

In the Computer Management window, click on the Groups folder to open it and select Action > New Group. Alternatively, use the right-click context menu.

The New Group dialog box appears.

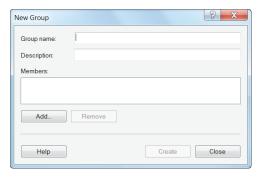

In the Group name field, enter the name of one of the groups specified in Table 8 on page 169 (TDS\_Administrator, TDS\_User, TDS\_Tech, TDS\_Guest). You can also enter a description in the Description field.


The group does not need any special operating-system level privileges.

- 3. Click Create to create and save the new group.
- Repeat this task for the remaining groups specified in Table 8 on page 169.

#### To add a user to a group on a local computer

In the New Group dialog box, click Add. Alternatively, double-click an existing group in the Groups folder to open its Properties dialog box, and click Add.

You see the Select Users dialog box.

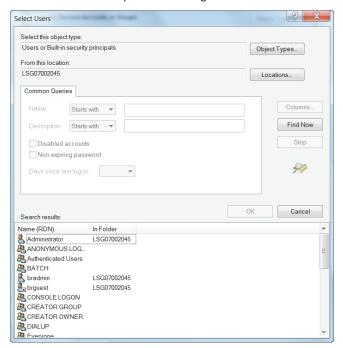

2. Click Advanced to expand the dialog box.

- In the expanded dialog box, click Find Now to populate the bottom field with all the users on the local computer.
- Click a user name in the list to select it, or press the Ctrl key and click multiple users to select them.
- 5. When you have selected all the users to add to the group, click OK, and click OK again to close the Select Users dialog box.
- 6. Click Create to close the New Group dialog box and create the group.
  - Alternatively, you can click OK to close the existing group's Properties dialog box and accept the changes.

## Configuring Users and Groups on a Network Domain

Note: The network administrator must know how users and groups are set up using the Windows server software at your site. The following example is used to illustrate the choices.

#### To locate the users and groups on a Windows server

Open Administrative Tools and select Active Directory.




Note that in the Active Directory window, the Users folder lists groups as well.

#### To create a new user on a Windows server

With the Users folder open, select Action > New User. Alternatively, use the right-click context menu.

You see the New User dialog box.

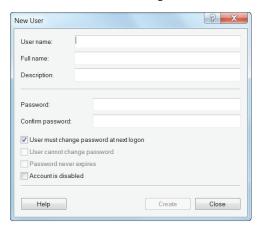

- 2. Fill out all the fields:
- **User name** The user name must be unique.
- **Full name** The full name must be unique.

Bio-Rad recommends using the user's actual full name, as this name will be shown in the document log and all the log reports. This is a requirement of 21 CFR 11.50a.

**Description** — This field must also be filled out.

Bio-Rad recommends that you enter the user's title as the description.

**Password** — Enter and confirm a password for the user.


Tip: Select the User must change password at next logon checkbox. This prevents the Windows system administrator from knowing the users' passwords.

**Note:** If you select the User must change password at next logon checkbox, the user must actually log on to Windows and change their password before using Image Lab Security Edition. Otherwise, Security Edition will not recognize the user.

### To create a new group on a Windows server

With the Users folder open, select Action > New Group. Alternatively, you can use the right-click menu.

You see the New Object - Group dialog box.

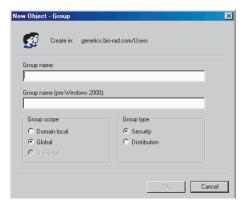

2. In the Group name field, enter one of the group names specified in Table 8 on page 169 (TDS\_Administrator, TDS\_User, TDS\_Tech, TDS\_Guest).


Type the name exactly as specified. You can also enter a description for the group in the Description field.

**Note:** The group does not need to have any special operating-system level privileges.

#### To add a user to a group on a Windows server

In the New Group dialog box, click Add. Alternatively, double-click an existing group in the User Manager folder to open its Properties dialog box and click Add.

You see the Select Users dialog box.

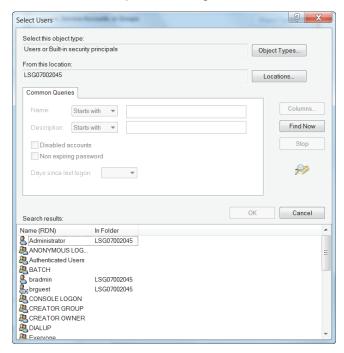

2. Click Advanced to expand the dialog box.

- In the expanded dialog box, click Find Now to populate the bottom field with all the users.
- Click a user name in the list to select it, or press the Ctrl key and click multiple users to select them.
- When you have selected all the users to add to the group, click OK, and click OK again to close the Select Users dialog box.
- 6. Click Create to close the New Group dialog box and create the group, or click OK to close the existing group's Properties dialog box and accept the changes.


## **Password Security**

21 CFR 11.300 (b) requires that passwords be "periodically checked, recalled, or revised." Password policies are therefore recommended, although the password duration and rules are up to the system administrator and the organization. For instance, the exact duration between password changes is flexible.


#### To set password policies on a local computer

- 1. Open the Control Panel and select Administrative Tools > Local Security Policy.
- 2. In the left pane of the Local Security Policy window, expand Account Policies and then select Password Policy.

You see the password policies listed in the right pane of the window.

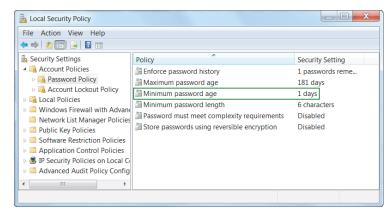

- 3. To change password policy settings:
  - Right-click the policy and select Properties to open its properties dialog box.

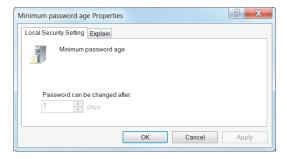

- Modify the default setting to meet your company policy.
- C. Click Apply.
- Click OK to close the Properties dialog box.
- Close the Local Security Policy window.

### To set password policies in Active Directory

- Open the Control Panel and select Administrative Tools > Domain Controller Security Policy.
- In the left pane of the Domain Controller Security Policy window, expand Security Settings > Account Policies and select Password Policy.
- 3. To change password policy settings:
  - Right-click the policy and select Properties to open its properties dialog box.
  - Modify the default setting to meet your company policy.
  - Click Apply. C.
  - Click OK to close the Properties dialog box.
- Close the Domain Controller Security Policy window.

## **Password Policy Setting Examples**

The following examples are only suggestions. Your organization needs to establish its own password policy.

- Enforce password history: 12 passwords remembered
- Minimum password age: 5 days
- Maximum password age: 30 days
- Minimum password length: 8 characters
- Password must meet complexity requirements: Enabled
- Store passwords using reverse encryption: Enabled

## **Account Lockout Policy Setting Examples**

- Account lockout duration: 0 (The account is locked out until the administrator unlocks it.)
- Account lockout threshold: 3 logon attempts
- Reset account lockout counter after: 30 minutes

## **Auditing Windows Event Logs**

Some global auditing information is stored in the Windows Event logs. It is a requirement of 21 CFR Part 11 that these logs be archived. However, by default, Windows systems automatically remove these data without warning.

Note: It is therefore critical that the event log is reconfigured to generate and preserve all necessary log data. Regular manual intervention is also required to preserve these data.

### To open the Event Properties Log

- Open Administrative Tools and click Event Viewer.
- 2. Right-click on each log and select Properties.

- In the section When maximum event log size is reached: select Do not overwrite events (Clear logs manually).
- 4. Increase the maximum size of the event log to cover any possible messages. The smaller the maximum size of the event log, the more often the system administrator must manually view, archive, and clear the system log.

Auditing information generated by the operating system is recorded in the Security Log. Logon failures in Image Lab Security Edition are recorded in this log.

During the review process, the log should be examined for attempted breaches of security, such as a series of failed logon attempts. To avoid the risk of losing data, the size should be very large and this inspection/archive process should occur daily. The Audit Policy should be set as follows:

- Audit account logon events Failure should be checked at a minimum
- Audit account management both Success and Failure should be checked
- Audit logon events Failure should be checked at a minimum
- Audit policy change both Success and Failure should be checked

## **Miscellaneous Security Measures**

Bio-Rad recommends that you use the built-in protections that Windows Server offers in order to protect the computer while the user is absent.

Note: Microsoft is continually updating its operating systems in response to security issues. It is critical to keep all components of the Windows operating system, especially any domain controllers, up to date.

## D Using Bio-Rad's **Stain-Free Technology**

Bio-Rad's stain-free gels eliminate the time-consuming staining and destaining steps required by other protein detection methods. Stain-free gels include unique trihalo compounds that allow rapid fluorescent detection of proteins with the Gel Doc™ EZ imager without staining.

When using Image Lab™ software version 5.1 or higher, the Gel Doc EZ imager is stain-free enabled to image these gels:

- Criterion™ TGX Stain-Free™ precast gels
- Criterion Stain Free™ precast gels
- Mini-PROTEAN® TGX Stain-Free™ precast gels
- TGX Stain-Free™ FastCast™ acrylamide solutions for handcast gels


When trihalo compounds in the gels encounter tryptophan residues, a UV light-induced reaction produces fluorescence, which can be easily detected by the Gel Doc EZ imager within gels or on low fluorescence polyvinyl difluoride (PVDF) membranes. Activation of the trihalo compounds in the gels adds 58 Da moieties to available tryptophan residues and is required for protein visualization. Proteins that do not contain tryptophan residues cannot be detected using this technology. The sensitivity of stain-free gels is comparable to staining with Coomassie Brilliant Blue for proteins with a tryptophan content >1.5%; sensitivity superior to Coomassie staining is possible for proteins with a tryptophan content >3%.

The benefits of stain-free technology include:

- Elimination of staining and destaining steps for faster time to results
- No background variability within a gel or between gels (as is often seen with standard Coomassie staining)
- Elimination of the need for acetic acid and methanol in staining and destaining, which reduces organic waste
- Visualization of transferred or blotted proteins on low fluorescence PVDF membranes

### Stain-Free Workflow

For detailed information about the Activate/image gels step, refer to Acquiring Images on page 69. For all other workflow steps, refer to the Criterion™ Precast Gels Instruction Manual and Application Guide (bulletin #4110001) or to the Mini-PROTEAN® Precast Gels Instruction Manual and Application Guide (bulletin #1658100).

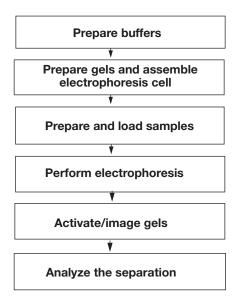

## **Electrophoresis with Stain-Free Gels**

Stain-free gels are made and packaged without sodium dodecyl sulfate (SDS), allowing them to be used for both SDS and native polyacrylamide gel electrophoresis (PAGE) applications.

#### To perform electrophoresis with stain-free gels

- 1. Prepare the sample and running buffers.
- 2. Set up the electrophoresis cell.
- Perform the run.

## **Imaging Gels**

Use unstained standards with stain-free gels, as some prestained standards are not compatible with stain-free technology. To monitor electrophoresis, use a 1:1 mixture of unstained and prestained standards.

Setting up a protocol for stain-free gels is very similar to setting up protocols for other applications. Follow the instructions in Setting Up a Custom Protocol on page 72. Choose one of the following activation times based on your sample and the purpose of your experiment:

- **Gels used in blotting** use 1 min activation for optimal results when performing western blotting followed by immunodetection.
- **Good sensitivity** use 2.5 min activation when samples are abundant and when a fully optimized signal-to-noise ratio is not necessary.
- **Best sensitivity** use 5.0 min activation for detection of proteins that are in low concentration and for the best quantification of the maximum number of bands. Because the reaction is near completion after 5 min, this method offers an optimal signal-to-noise ratio.

Note: If the gel has already been activated for 2.5 min, activating it for another 2.5 min might improve it. But activating an image for more than 5 min will not.

## **Imaging Blots**

To blot stain-free gels, use standard blotting procedures as described in the instruction manual you are using. Use only PVDF membranes with low background fluorescence, as membranes other than low fluorescence PVDF can result in high background or low sensitivity with the Gel Doc EZ imager.

To assess transfer efficiency, be sure to activate and visualize the gel using the imager before transfer.

## **Regression Calculation** Methods

Each regression method calculates a standard curve. Some of the methods provide the formula for the standard curve. In this case, the molecular weight can be calculated by:

x = relative front of the band of interest

y = molecular weight of the band of interest

**Linear (semilog):** The linear equation is y = a + bx, where a is the intercept and b is the slope of the line.

Note: The linear equation is calculated on the log of the molecular weight values.

The R<sup>2</sup> value can be used to determine the overall quality of the linear fit. A linear regression with an R2 value of >0.99 is considered a very good fit. The primary advantage of this method is that it is extremely simple. The primary disadvantage is that it will deliver incorrect results if the data are not very linear.

Point-to-point (semilog): No single equation is available for the point-to-point method. The slope of each segment of the curve between data points is calculated independently.

Note: The log of the molecular weight values is used to calculate the slope for each segment of the curve.

Logistic: The logistic-4PL equation is

$$y = d + \frac{a - d}{1 + \left(\frac{x}{c}\right)^b}$$

where:

x = mobility

y = molecular weight

a = estimated molecular weight at infinity

b = slope of the tangent at midpoint

c = midpoint

d = estimated molecular weight at zero mobility

Since the curve generated by the logistic-4PL regression method represents a perfectly shaped S, it might not fit the data very well in all cases.

Cubic spline: Cubic spline curves are smooth curves that go through every data point. The model is a cubic polynomial on each interval between data points. In some cases, a spline curve can work well as a standard curve for interpolation. However, because the curve is calculated individually for every pair of points, it does not correspond to any single equation.

## Mitsubishi P93/P95 **Thermal Printer**

## **Setting up a Thermal Printer on Windows**

The printer driver is on the Image Lab™ software installation CD in the Misc folder.

#### To set up a thermal printer on a Windows system

- 1. Install the printer driver.
- Open the printer section in Control Panel.
- Click the thermal printer icon and select Printing Preferences.



- Configure the correct paper size. Select 1280 x 1280 from the dropdown list.
- Click OK to apply your changes.

## Setting up a Thermal Printer on a Mac

The printer driver can be found on the Image Lab software installation CD in the Misc directory.

#### To set up a thermal printer on a Mac system

- 1. Install the printer driver.
- 2. Connect the printer to the computer.

#### To configure the correct paper size

- 1. Start Image Lab software.
- 2. Select File > Page Setup.



- In the Settings dropdown list, select Page Attributes.
- 4. In the Format For dropdown list, select the Mitsubishi printer.
- 5. In the Paper Size dropdown list, select 1280 x 1280.
- 6. In the Settings dropdown list, select Save as Default.
- 7. Click OK to save the settings.

# **Glossary**

Aspect ratio	The ratio of the width to the height of an image.
CCD	(Charge-coupled device) A light-sensitive silicon chip used as a photodetector in ChemiDoc™ MP and ChemiDoc XRS+ camera systems.
Colormaps	Different color representations of a gel image.
Electrophoresis	A technique for separating molecules based on the differential movement of charged particles through a matrix when subjected to an electric field.
Example precision	The number of decimal places chosen for displaying a measurement.
Flat fielding	An average intensity computation that compensates for nonuniformities generated by an instrument.
Histogram	A graphed representation of the brightness, or gray value, of an image.
Native charge density	The inherent electrical charge of a protein without the addition of SDS.
pl	Isoelectric point; the pH at which a protein molecule carries no net charge.
Rf	Relative front value of the band. In Image Lab™ software, Rf has a value between 0 and 1 and indicates the relative movement of the band from top to bottom.
Quantitative imaging	Determines the quantity of a protein's components through analysis of the pixel values in a digital image of the sample.
UV-B	The range of ultraviolet light used by the system.
UV transilluminator	The part of the imager that transmits UV light through a sample.





Bio-Rad Laboratories, Inc.

Life Science Group Web site www.bio-rad.com USA 800 424 6723 Australia 61 2 9914 2800 Austria 01 877 89 01 Belgium 09 385 55 11 Brazil 55 11 3065 7550 Canada 905 364 3435 China 86 21 6169 8500 Czech Republic 420 241 430 532 Denmark 44 52 10 00 Finland 09 804 22 00 France 01 47 95 69 65 Germany 089 31 884 0 Greece 30 210 9532 220 Hong Kong 852 2789 3300 Hungary 36 1 459 6100 India 91 124 4029300 Israel 03 963 6050 Italy 35 02 216091 Japan 81 3 6361 7000 Korea 82 2 3473 4460 Mexico 52 555 488 7670 The Netherlands 0318 540666 New Zealand 64 9 415 2280 Norway 23 38 41 30 Poland 48 22 331 99 99 Portugal 351 21 472 7700 Russia 7 495 721 14 04 Singapore 65 6415 3188 South Africa 27 861 246 723 Spain 34 91 590 5200 Sweden 08 555 12700 Switzerland 026674 55 05 Taiwan 886 2 2578 7189 Thailand 1800 88 22 88 United Kingdom 020 8328 2000

10019683 Rev F US/EG 13-2298 1213 Sig 1213